






Determine number of fluorines attached to each carbon by ^{13}C NMR spectroscopy!

The number of fluorine attached to each carbon is useful in structure analysis of fluorinated compounds. ^{19}F signals are sometimes observed in a very wide chemical shift range. In that case, It is difficult to uniformly excite all ^{19}F signals, and hence APT(Attached Proton Test) experiment is more useful than DEPT.

On the other hand, APT is less sensitive than DEPT, so we need to set a higher number of scans. The figures below show ^{19}F , ^{13}C and APT spectra of 5% ethyl 1,1,2,3,3,3-hexafluoropropyl ether in $\text{DMSO-}d_6$. This sample has ^{19}F signals spread over a frequency range of over 50 kHz, and so uniform excitation is very challenging.

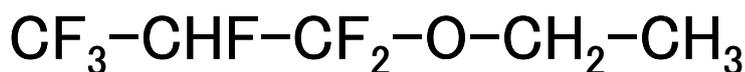

- In the ^{13}C - ^{19}F APT spectra, ^{13}C signals of CF_3 and CF have positive signals, while CF_2 , C and solvent signals have negative signals.
- By the application of ^1H decoupling, we can increase sensitivity!
- ROYALPROBE HFX can perform these ^1H , ^{19}F , ^{13}C triple-resonance measurements even with a standard 2-channel console!

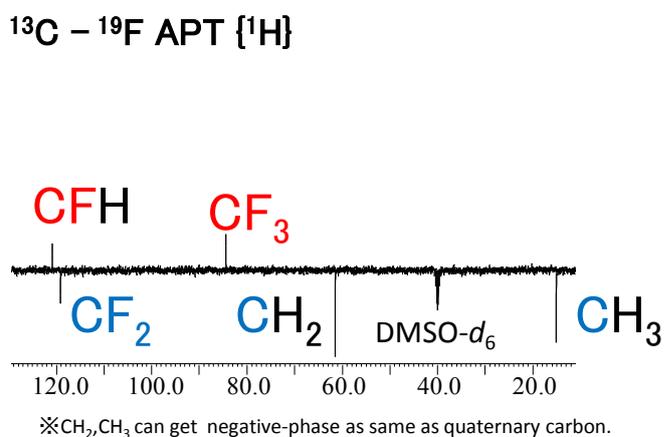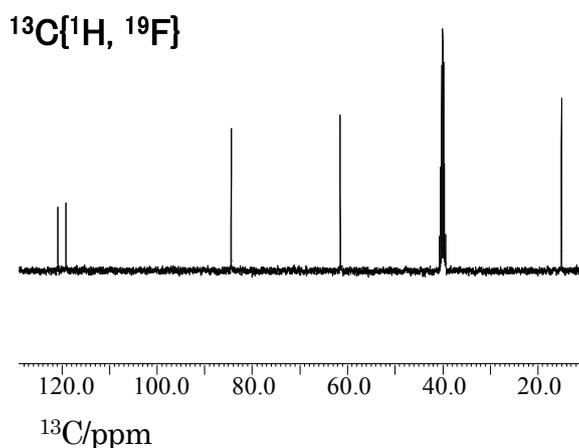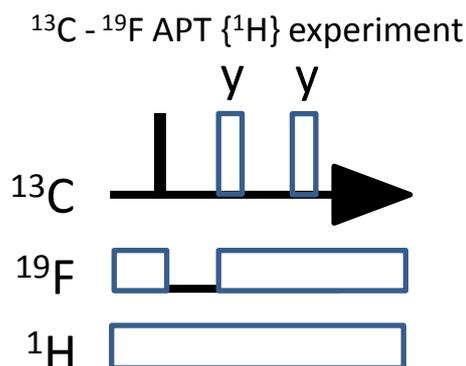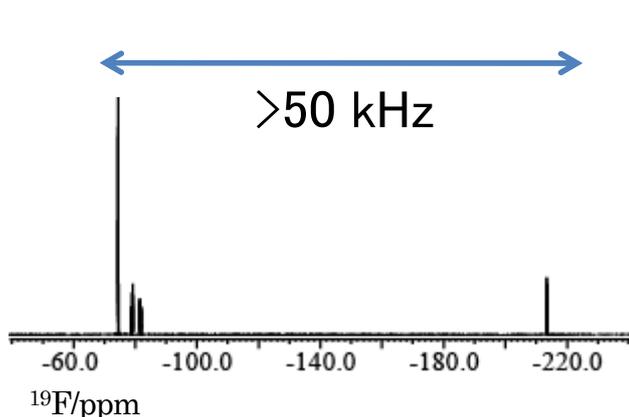

$^{13}\text{C}\{^1\text{H}, ^{19}\text{F}\}$ and ^{13}C - ^{19}F APT $\{^1\text{H}\}$ spectra, 32 scans

console : JNM-ECZ400S, ROYALPROBE HFX

Copyright © 2018 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.

