

Delta Tips

NMDT_0035


Printing: How to Remove the JEOL Logo

NMR data processing software

Delta
NMR Software
v5.0




As a default, the JEOL logo is printed whenever you print a spectrum to a printer, PDF file or clipboard. If you wish not to print the logo, follow the procedure described in this issue.

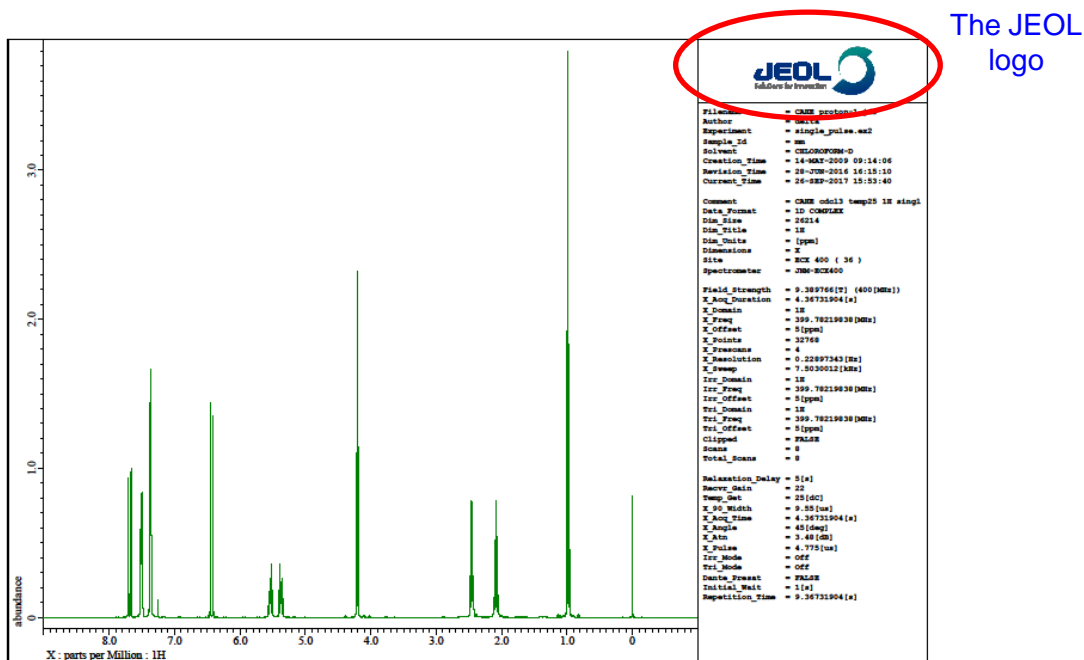


① Select **Options – Preferences** in the Delta window.

The Delta window


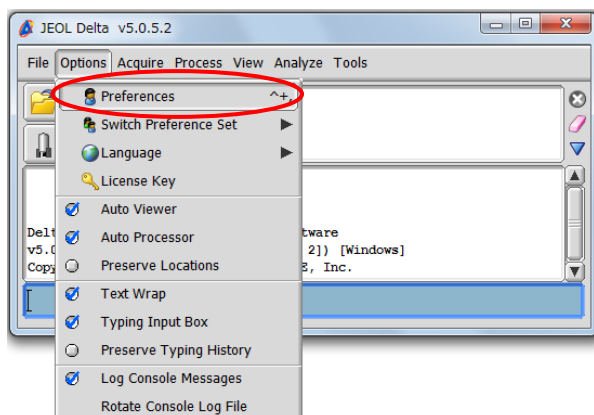


The Preferences window

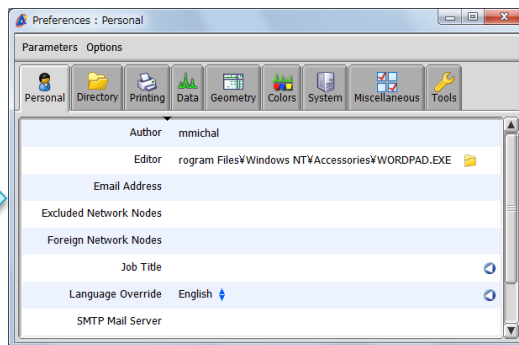

Delta Tips

- ② Select the **Printing** tab.
- ③ Turn off the **Plot JEOL Logo** option and close the **Preferences** window.

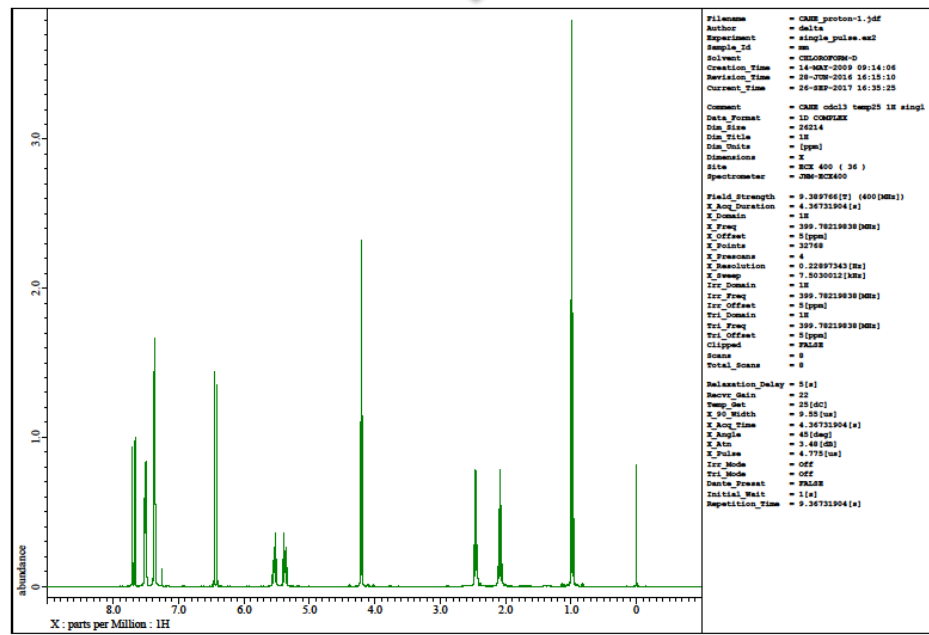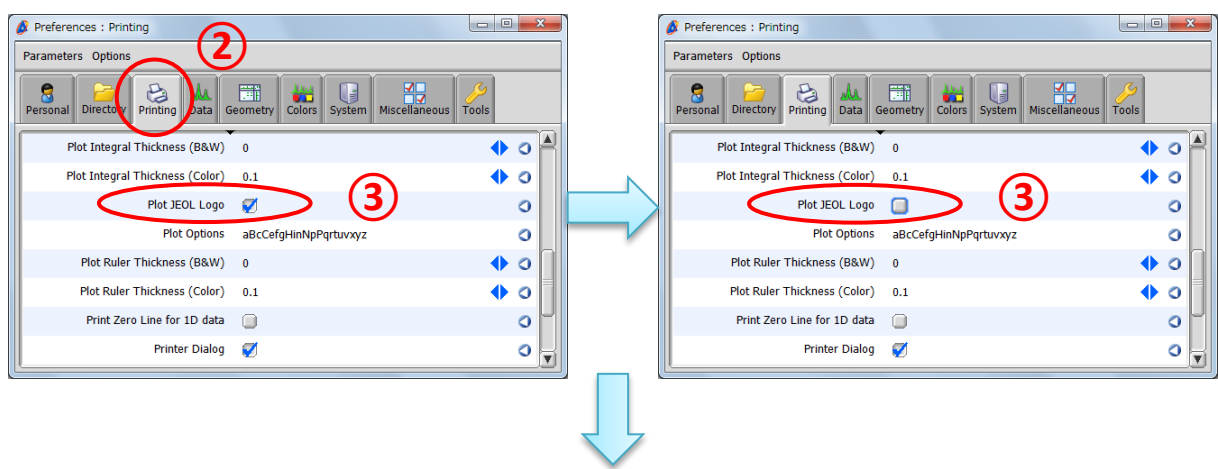

! The modification of **Preferences** is permanent.

★ If you wish to reset any parameter of **Preferences** to its default value, click the  button.

