

Delta Tips

NMDT_0043

Data Slate (Part 6)

NMR data processing software


Delta
NMR Software
v5.0



< How to align multiple spectra variously >

Data Slate can display multiple spectra in various ways. 1D and/or 2D spectra can be aligned horizontally, vertically and in boxes.


★ Refer to NMDT_0025 for the display function of multiple spectra in one Data Slate window.

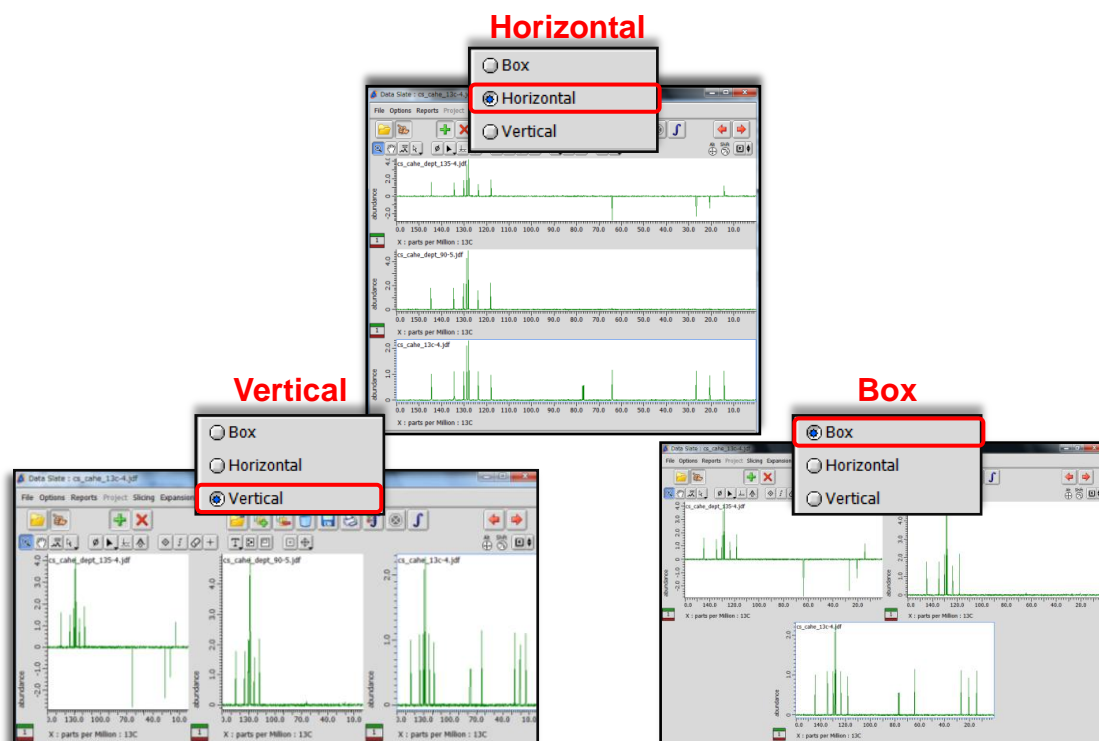

Select **Options – Box (Horizontal or Vertical)** to change the alignment of the multiple spectra.

In the example below, three spectra are displayed in the **Box** mode.

Options menu

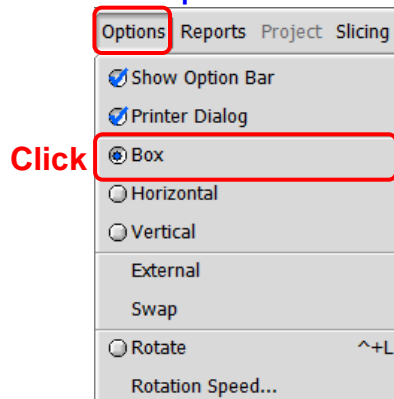

The Box alignment of multiple spectra

