

Simplifying ^1H -NMR Spectra in NMR Metabolic Profiling —Applying Projections of ^1H homonuclear 2D J -resolved spectra—

Product used : Nuclear Magnetic Resonance (NMR)

Spectral overlap often hampers subsequent analyses (ex. multivariate analyses) in ^1H -NMR spectra of molecular mixtures such as in NMR metabolic profiling (NMR-MP). One of the most straightforward solution is using more higher field NMR system. However it is not always easy to address. In this note, the method simplifying ^1H -NMR spectra of mixture is introduced by use of projections of ^1H homonuclear 2D J -resolved spectra.

Simplifying ^1H -NMR spectrum by applying projections of ^1H homonuclear 2D J -resolved spectrum [1]

In the 2D J -resolved spectroscopy experiment, an evolution of J -coupling is encoded while that of chemical shift is refocused in the indirect evolution period. Therefore the proton multiplets appear at a 45 degree angle in the resultant 2D spectrum as shown in **Figure 1(a)**. Then this 2D spectrum is tilted by 45 degrees to remove J -couplings from x-axis. After tilting, as shown in **Figure 1(b)**, the 2D spectrum contains chemical shift and J -coupling alone in x- and y-axis, respectively. The projection of the x-axis is a pseudo pure shift spectrum (**Figure 1(c)**: upper). You can see how spectrum goes is simplified relative to a standard single pulse experiment.

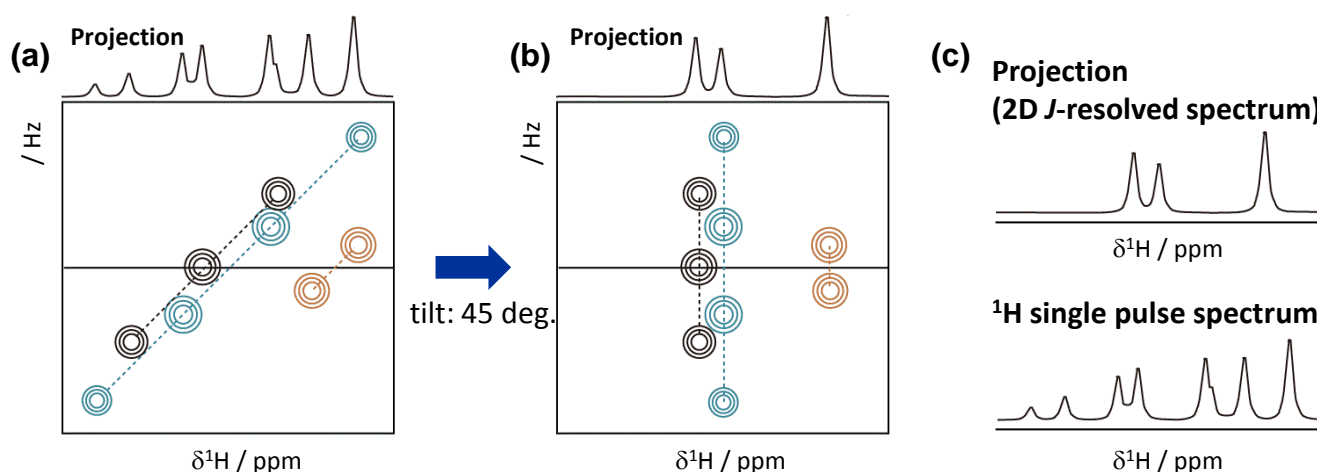


Figure 1 Simplifying ^1H -NMR spectrum by applying projections of ^1H homonuclear 2D J -resolved spectrum. (a) The proton multiplets appear at a 45 degree angle in original 2D spectrum. (b) After tilting, the 2D spectrum contains chemical shift and J -coupling in x- and y-axis, respectively. (c) The projection of (b) is more simple than the conventional ^1H single pulse spectrum.

Suppressing artifacts from strong couplings in J -resolved spectroscopy

We can record simple spectra using projections of 2D J -resolved spectra even though standard magnetic field NMR system such as 400 MHz system is used. However there is a drawback in J -resolved spectroscopy that artifacts from strong coupling can occur, especially when a lower magnetic field system is used.

As it was introduced before in our application note "Suppressing artifacts from strong couplings in 2D J -resolved spectroscopy" (NM050018), these artifacts can be suppressed when double spin-echo type of 2D J -resolved spectroscopy [2] is used.

A projection of 2D J-resolved spectrum of polar metabolites in broccoli sprout.

There are often severe spectral overlap in ^1H -NMR spectra when plants are used as samples for NMR-MP because they produce various kind of metabolites including plant specialized metabolites. Here, we demonstrate that using projections of 2D J-resolved spectra is a powerful method in actual NMR-MP studies. Polar metabolites in broccoli sprouts were extracted and used as a NMR sample. All spectra were recorded with standard NMR system (JNM-ECZ400S spectrometer with a ROYALPROBE). The double spin-echo method was used as a 2D J-resolved spectroscopy experiment.

^1H spectra of polar metabolites in broccoli sprout are shown in **Figure 2**. **Figure 2(a)** and **(b)** show a projection of 2D J-resolved spectrum and ^1H single pulse spectrum, respectively. As many metabolites were present in the sample, there is severe spectral overlap. Using a projection of 2D J-resolved spectrum, we were able to record far simpler spectra than the results of a typical ^1H single pulse spectrum. For example, the insets in the spectra below highlight four different methyl protons in valine and isoleucine that were clearly resolved in the projection of 2D J-resolved spectrum **(a)** whereas they were ambiguous in ^1H single pulse **(b)**.

2D J-resolved spectroscopy is also useful to assign metabolites [3]. Please see detail of an actual NMR-MP study using projections of 2D J-resolved spectra in our application note "Unsupervised Multivariate Analyses for NMR-based Metabolomics" (NM170013) [4].

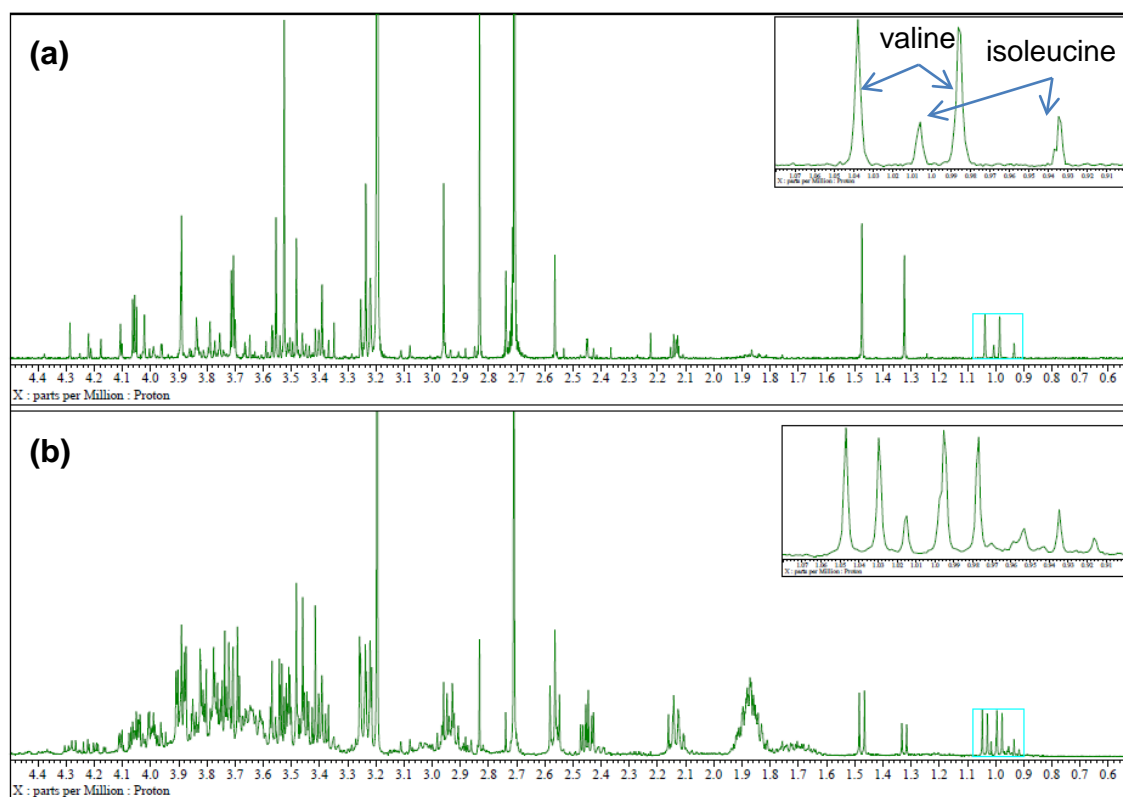


Figure 2 A comparison between a projection of 2D J-resolved spectrum and conventional ^1H single pulse spectrum of polar metabolites in broccoli sprout. (a) A projection of 2D J-resolved spectrum. The double spin-echo method was used. (b) ^1H single pulse spectrum.

References

- [1] Ludwig; C., Viant; M., *Phytochem. Anal.*, **21** (2010) 22-32.
- [2] Thrippleton; M. J., Edden; R. A. E., Keeler; J., *J. Magn. Reson.*, **174** (2005) 97.
- [3] Kikuchi; J. et al., *Anal. Chem.*, **88** (2016) 659.
- [4] Application note "Unsupervised Multivariate Analyses for NMR-based Metabolomics" (NM170013).

Copyright © 2017 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.



JEOL Ltd.

3-1-2 Musashino Akishima Tokyo 196-8558 Japan Sales Division Tel. +81-3-6262-3560 Fax. +81-3-6262-3577
www.jeol.com ISO 9001 • ISO 14001 Certified

• **AUSTRALIA & NEW ZEALAND** /JEOL(AUSTRALASIA) Pty.Ltd. Suite 1, L2 18 Aquatic Drive - Frenchs Forest NSW 2086 Australia • **BELGIUM** /JEOL (EUROPE) B.V. Planet II, Gebouw B Leuvensesteenweg 542, B-1830 Zaventem Belgium • **BRAZIL** /JEOL Brasil Instrumentos Científicos Ltda. Av. Jabaquara, 2958 5º andar conjunto 52 ; 04046-500 Sao Paulo, SP Brazil • **CANADA** /JEOL CANADA, INC. 3275 1ere Rue, Local #8 St-Hubert, QC J3Y-8Y6, Canada • **CHINA** /JEOL(BEIJING) CO., LTD. Zhongkeziyuan Building South Tower 2F, Zhongguancun Nanshanjie Street No. 6, Haidian District, Beijing, P.R.China • **EGYPT** /JEOL SERVICE BUREAU 3rd Fl. Nile Center Bldg., Nawal Street, Dokki, (Cairo), Egypt • **FRANCE** /JEOL (EUROPE) SAS Espace Claude Monet, 1 Allée de Giverny 78290, Croissy-sur-Seine, France • **GERMANY** /JEOL (GERMANY) GmbH Gute Aenger 30 85356 Freising, Germany • **GREAT BRITAIN & IRELAND** /JEOL (U.K.) LTD. JEOL House, Silver Court, Watchmead, Welwyn Garden City, Herts AL7 1LT, U.K. • **ITALY** /JEOL (ITALIA) S.p.A. Palazzo Pacinotti - Milano 3 City, Via Ludovico il Moro, 6/A 20080 Basiglio(MI) Italy • **KOREA** /JEOL KOREA LTD. Dongwoo Bldg. 7F, 1443, Yangjae Daero, Gangdong-Gu, Seoul, 05355, Korea • **MALAYSIA** /JEOL(MALAYSIA) SDN.BHD. 508, Block A, Level 5, Kelana Business Center, 97, Jalan SS 7/2, Kelana Jaya, 47301 Petaling Jaya, Selangor, Malaysia • **MEXICO** /JEOL DE MEXICO S.A. DE C.V. Arkansas 11 Piso 2 Colonia Napolis Delegacion Benito Juarez, C.P. 03810 Mexico D.F., Mexico • **RUSSIA** /JEOL (RUS) LLC. Krasnoproletarskaya Street, 16, Bld. 2, 127473, Moscow, Russian Federation • **SCANDINAVIA** /SWEDEN JEOL (Nordic) AB Hammarbacken 6A, Box 716, 191 27 Sollentuna Sweden • **SINGAPORE** /JEOL ASIA PTE.LTD. 2 Corporation Road #01-12 Corporation Place Singapore 618494 • **TAIWAN** /JIE DONG CO., LTD. 7F, 112, Chung Hsiao East Road, Section 1, Taipei, Taiwan 10023 (R.O.C.) • **THE NETHERLANDS** /JEOL (EUROPE) B.V. Lieweg 4, NL-2153 PH Nieuw-Vennep, The Netherlands • **USA** /JEOL USA, INC. 11 Dearborn Road, Peabody, MA 01960, U.S.A.