

Delta Tips

NMDT_0006

User Authentication

NMR data processing software

Delta
NMR Software
v5.0



When connecting to a spectrometer, you can type the user name or select it from the list of users after pressing the  button. However, if the Delta user name is identical to the current Windows user name, there is a way to load the user name with a single click. In addition to this simple trick, if the Delta user name is identical to the Delta password, it can be copied from the Name input box to the Password input box.

- ① Click the  button in the **Authentication** window to load the user name.

Authentication

Please enter your login information

Name Account name required

Password

Connect Own Cancel

- ② Place the cursor inside the **Name** input box, press the right mouse button and drag & drop 'delta' to the **Password** input box.

Authentication

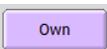
Please enter your login information

Name delta

Password

Connect Own Cancel

Drag & drop 'delta'

- ③ Click either the  or  button to connect to the spectrometer.

Authentication

Please enter your login information

Name delta

Password *****

Connect Own Cancel