

# Delta Tips

NMDT\_0011

## How to Open Periodic Table Tool


NMR data processing software

Delta  
NMR Software  
v5.0



The Periodic Table Tool displays NMR related information, such as natural isotopes, natural abundance, nuclear spin number, resonance frequency and relative sensitivity.

Select **Tools – Periodic Table** in the JEOL Delta, Data Slate or Processor window.

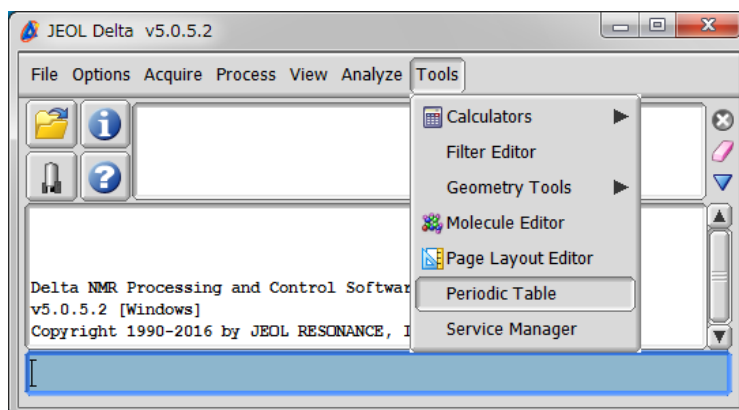

- ① Input  $^1\text{H}$  resonance frequency. The default value is 100 MHz.
- ② Select an element.

Legend

Isotope	Abundance	Spin	Frequency	Rel. 1H	Rel. 13C
$^{14}\text{N}$	99.63	+1	28.904	0.00101	5.69
$^{15}\text{N}$	0.37	-1/2	40.544	0.00104	0.0219