

Delta Tips

NMDT_0018

Curve Analysis Tool

NMR data processing software


Delta
NMR Software
v5.0




< How to plot relaxation or diffusion data of multiple peaks at once >

The procedure is demonstrated on T_1 relaxation data.

- Click the  button to activate the selection mode. Select the peaks of interest to plot their relaxation data.

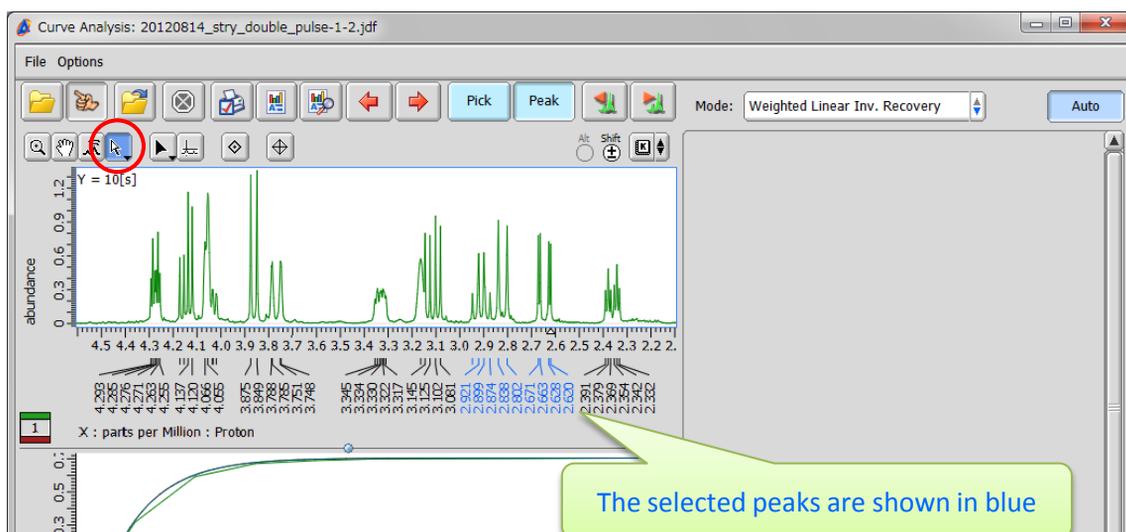


- Click the  button to plot the data. The **Print Options** window opens.

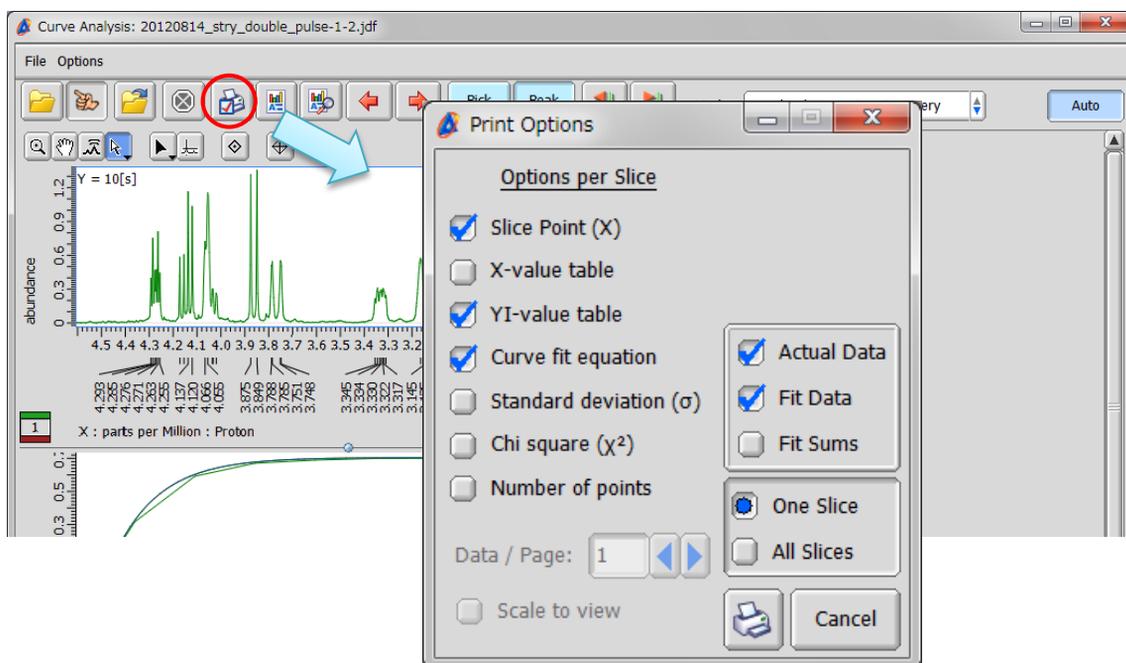


- ③ Select the 'All Slices' option in the Print Options window.
- ④ It is possible to print up to four plots per page. We will plot four plots per page.

The Print Options window

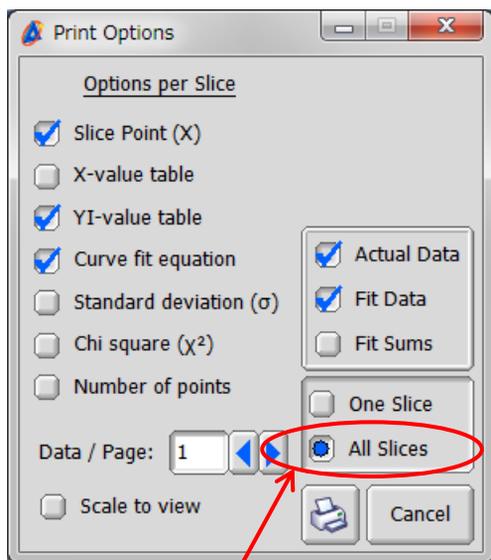

③ Print data for multiple peaks

The Print Options window

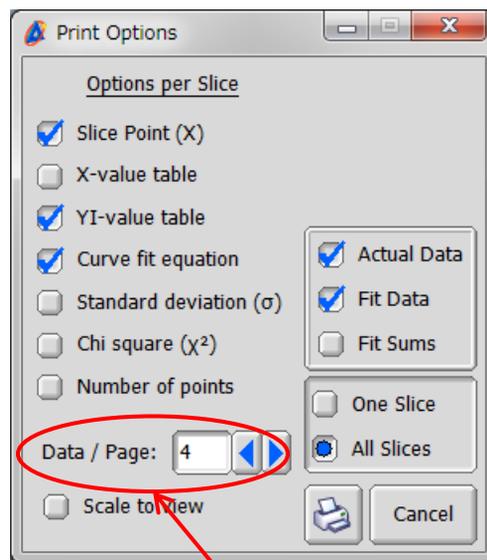

④ The number of plots per page

- ⑤ Click the  button in the **Print Options** window to plot the data. In the **Print** window (not shown), select a printer to plot on paper or JEOL PDF to print to a PDF file.

- ⑥ An example of plot is shown below.

