

Delta Tips

NMDT_0023

How to Export Spectrum to ASCII File

NMR data processing software


Delta
NMR Software
v5.0




Open the data file to export to the ASCII format with 1D Processor or Data Slate.

- ① Select **File – Save As (1D Processor)** or **File – Save – Save As (Data Slate)**. Alternatively, push the **Shift, Ctrl and S** keys simultaneously.


The 1D Processor window

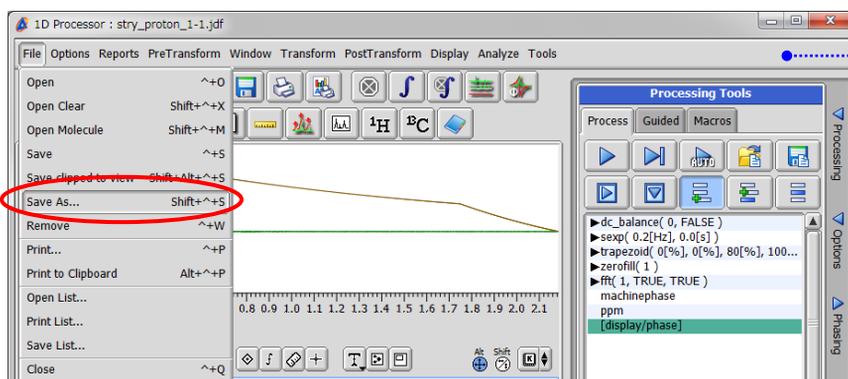

The Data Slate window

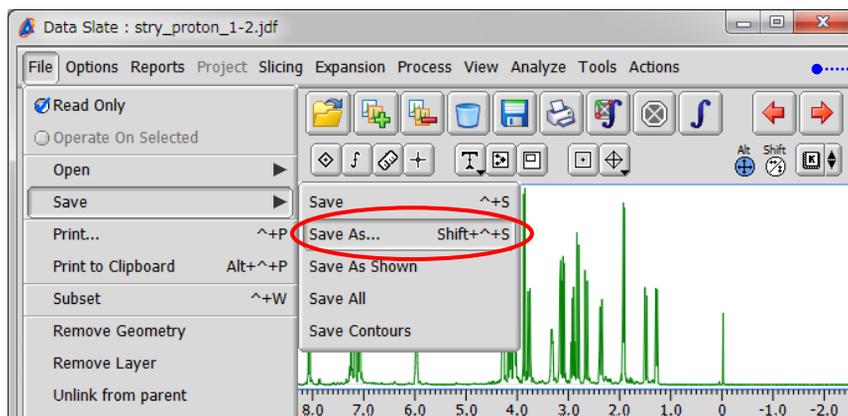


- ② Select **File – Export Format – JEOL : Generic ASCII** and click the  button.

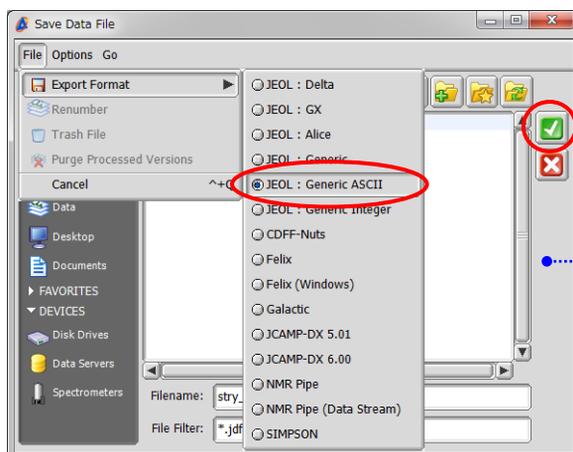

The Save Data File window



Delta Tips

- ③ The conversion process starts and an **Inform** window appears. After the process has finished, the window automatically closes and a report appears in the Delta window.

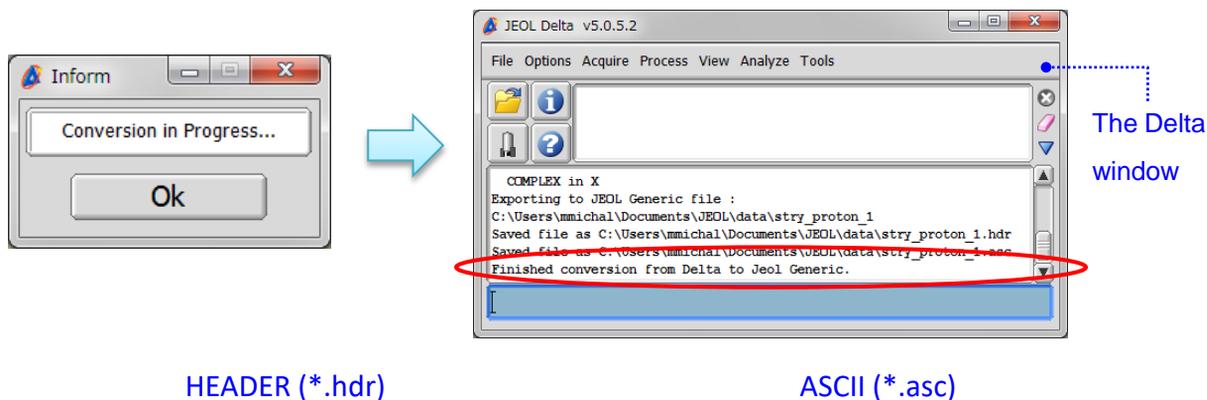

HEADER (*.hdr)

ASCII (*.asc)

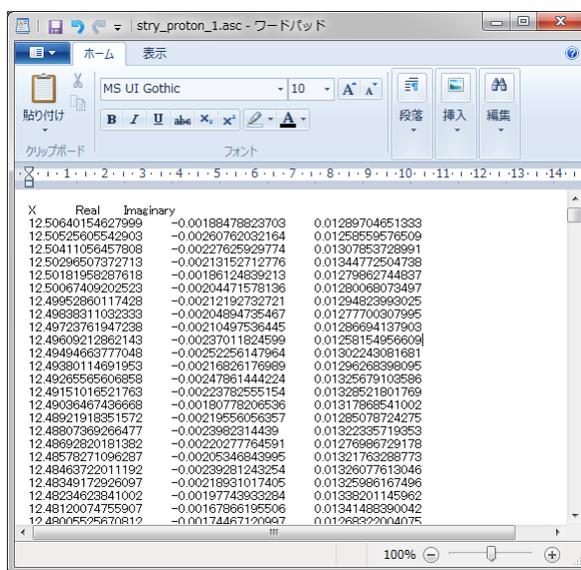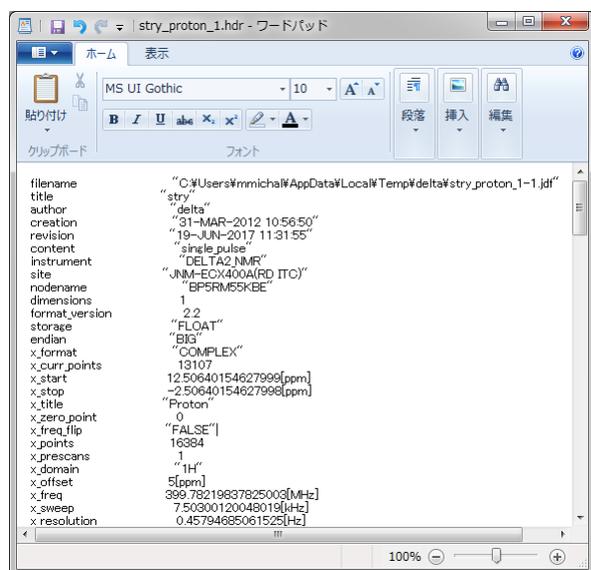

- ★ If you wish to export FID data into the ASCII format, use **1D Processor**. However, delete the processing list before following STEPS 1 – 3.

