

Delta Tips

NMDT_0024

Picture-in-Picture (PIP) Function


NMR data processing software


Delta
NMR Software
v5.0



< How to hide the box indicating the range in 1D spectrum >

The PIP function allows you to display a spectrum while zooming up range/signal of interest.

To create a PIP box click the  icon.

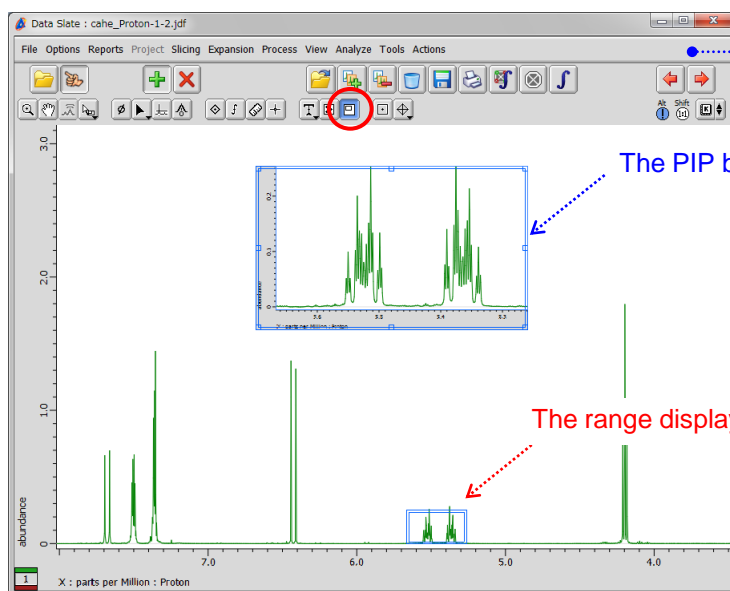

If you wish to hide the blue box indicating the range displayed in the PIP box, push and hold a right-mouse-button and unselect **Options – PiP Windows**. Alternatively, push the **Alt** and **W** keys simultaneously.

