

# Delta Tips

NMDT\_0027

## Combine Files Tool

NMR data processing software

Delta  
NMR Software  
v5.0





### < How to use Combine Files Tool >

When the Combine Files Tool is used, it is possible to combine multiple 1D data files and save them as one pseudo-2D data file. It is possible to combine spectra or FIDs (Free Induction Decay). We will combine five individual proton spectra in this example.

- ① Select **Process – Combine Files** to open the Combine Files tool.

The Data Slate window

The Combine Files window




- ② Click the  icon and then  to select a file in the fingering mode. Select all the spectra one by one.

The Combine Files window


# Delta Tips



③ Specify the number of files (spectra) to combine as shown in the figure. It is 5 in our example.

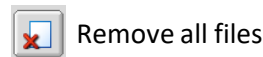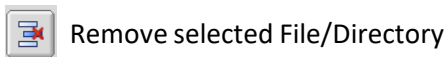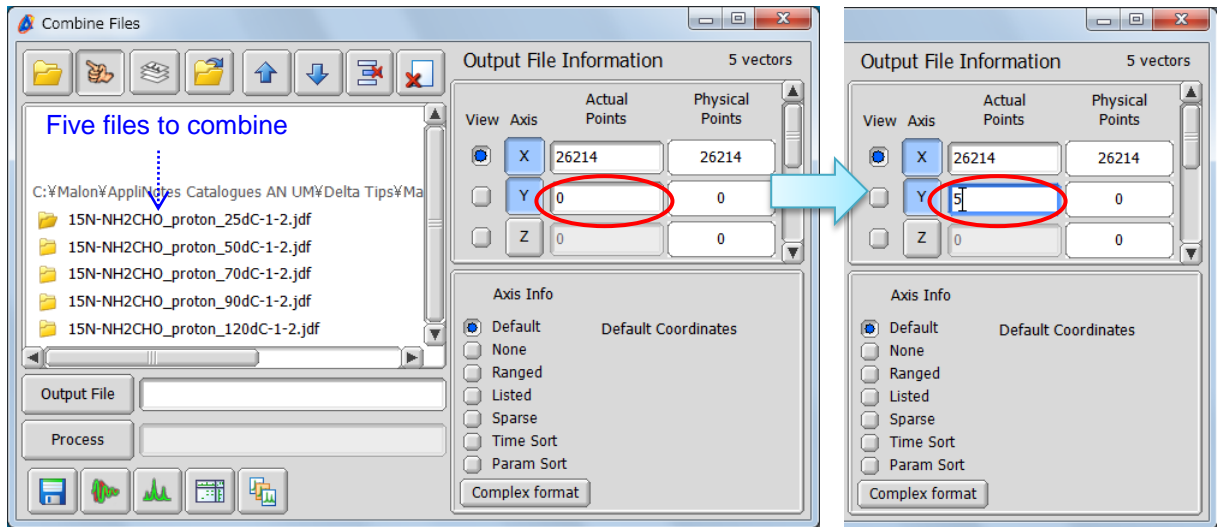


④ Type a filename into the Output File box.

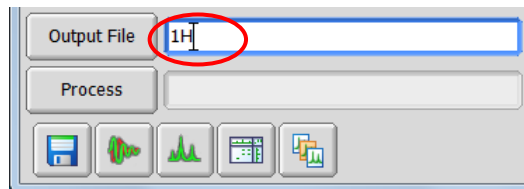

⑤ Click the  button to save the new file.

★ If you click the  button, a new window opens in which you may specify location.

You can click the  button and switch to the Stack Plot mode.

