

# Delta Tips

NMDT\_0034

## Precision of Chemical Shifts and Integrals

NMR data processing software

**Delta**  
NMR Software  
v5.0




< How to adjust the number of digits after the decimal point >


The precision (scale) of chemical shifts and integrals can be adjusted via **Preferences**.

Select **Options – Preferences** in the Delta window.

The Delta window

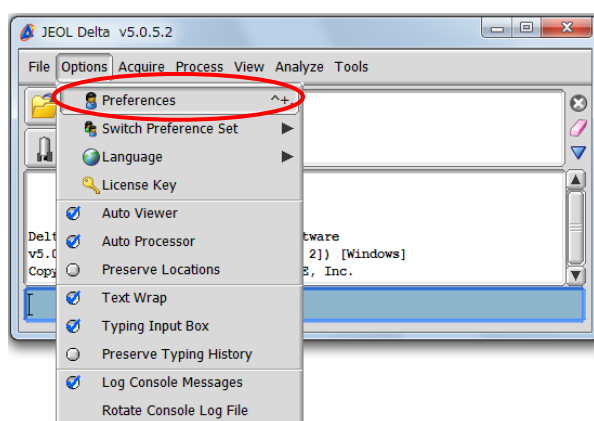

The Preferences window

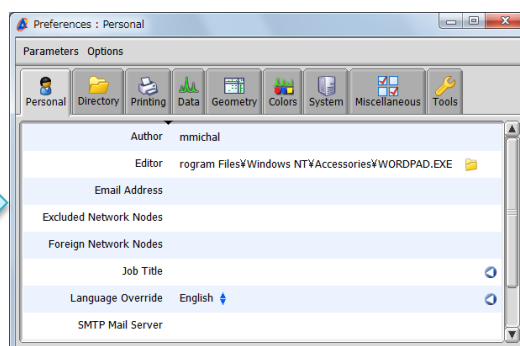


! The modification of **Preferences** is permanent.

## 1 Precision of Chemical Shifts


- ① Select the **Geometry** tab.
- ② Modify the **Peak Precision** value and close the **Preferences** window.

① Geometry tab

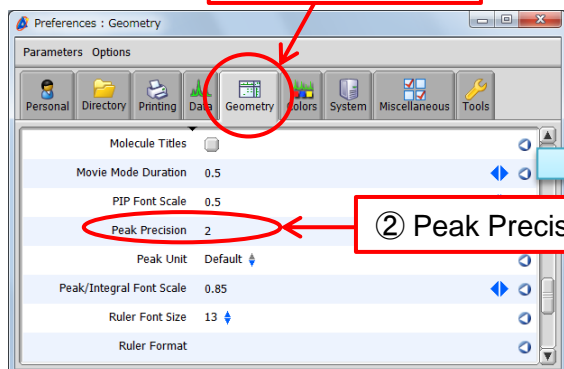


Two decimal places are shown

② Peak Precision

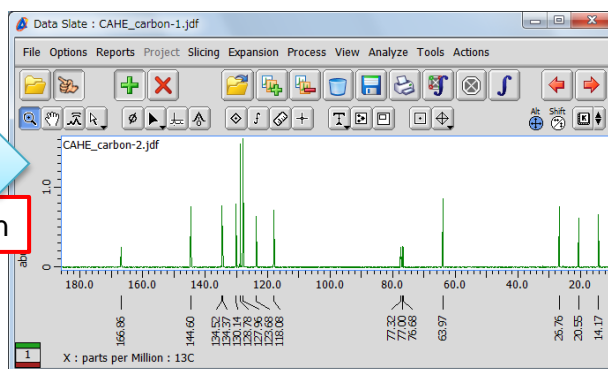

## 2 Precision of Integrals

- ① Select the **Geometry** tab.
- ② Modify the **Integral Format** value and close the **Preferences** window.

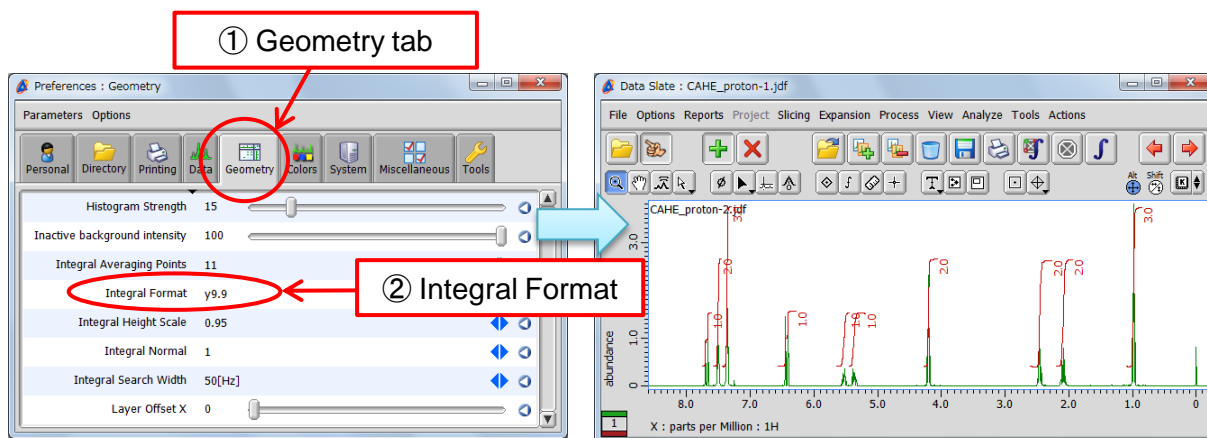

One decimal place is shown



y9.99 → Two digits after the decimal point  
y9.9 → One digit after the decimal point