

# Delta Tips

NMDT\_0041

## How to display/hide Y-axis ruler


NMR data processing software

Delta  
NMR Software  
v5.0



It is possible to instantly display or hide the Y-axis ruler in the Data Slate or Processor window.


Push the **Alt** and **Y** keys to activate/deactivate the Y-axis ruler display function.

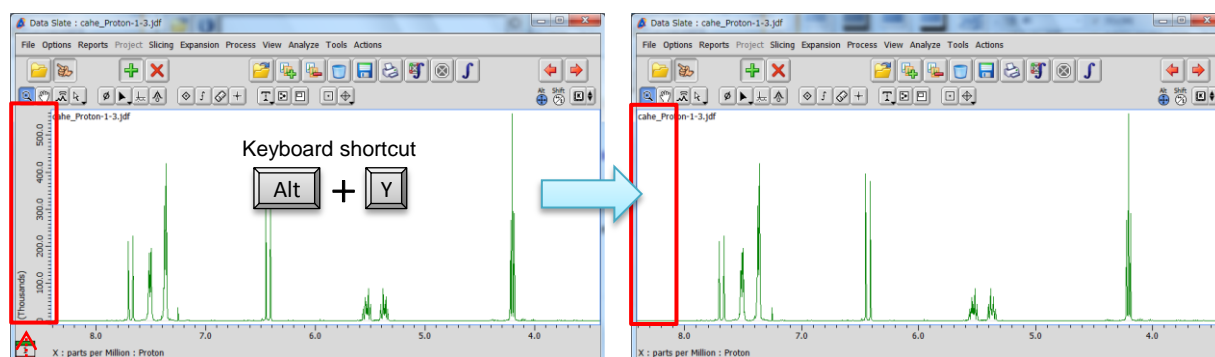

Y-axis ruler

Note that the Y-axis ruler has been hidden

★ It is also possible to display or hide Y-ruler from the context menu as follows:  
Push and hold the right mouse button to display the menu. Then select **Options – Rulers – Y Ruler** as shown below.

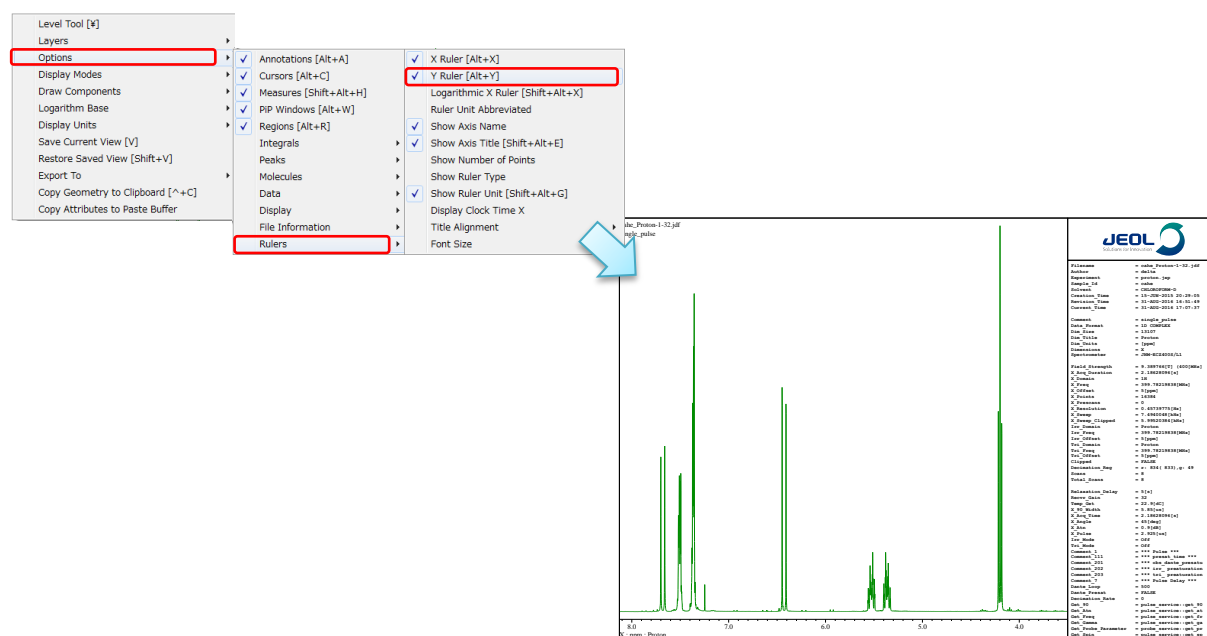

An example of <sup>1</sup>H NMR spectrum without Y-ruler