

Delta Tips

NMDT_0044

Data Slate (Part 7)

NMR data processing software

Delta
NMR Software
v5.0



< How to swap one spectrum for another one >

It is possible to exchange position of one spectrum for position of another spectrum.

★ Refer to NMDT_0025 and 0043 for the display functions of multiple spectra in one Data Slate window.

① Select **Options – Swap** to activate the Swap functions. The cursor has changed into **Finger** symbol .

② Select the spectra to swap with the cursor. In the example below, we have exchanged the position of spectrum **A** for the position of spectrum **C**.

