

Delta Tips

NMDT_0045

Data Slate (Part 8)

NMR data processing software


Delta
NMR Software
v5.0



<How to open a spectrum in another Data Slate window>

The Data Slate window has the function of opening a spectrum in another Data Slate.

- ★ Refer to NMDT_0025 for the display function of multiple spectra in one Data Slate window.

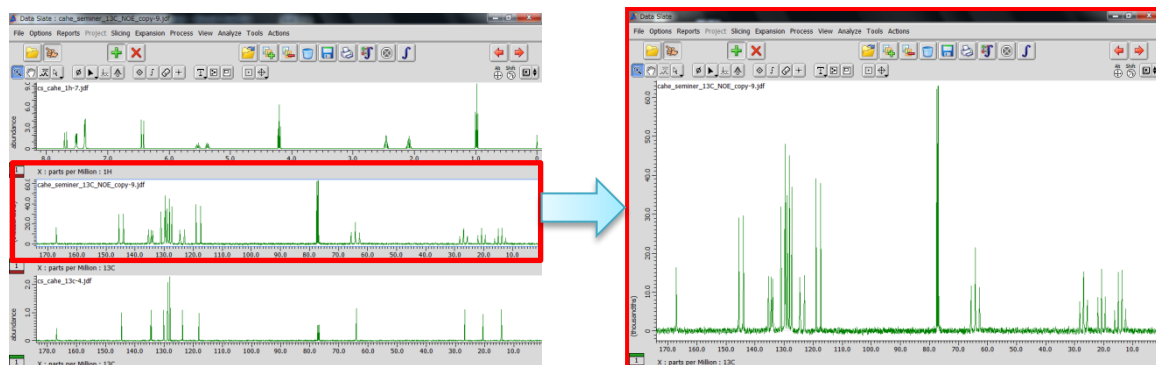

- ① Select the spectrum which you wish to open in a new Data Slate window.
- ② Select **Options – External** to open the selected spectrum in the new Data Slate window.

In the example below, we have used the middle spectrum (DEPT-135).

Options menu

The DEPT-135 spectrum has opened in the new window

