

Delta Tips

NMDT_0048

Peak Sequence Markers

NMR data processing software

Delta
NMR Software
v5.0



It is possible to display peak sequence markers on 2D data in the **2D Viewer** window.


You can active/deactivate the peak sequence markers function by pushing the



[Alt+M] keys.

To reproduce the example demonstrated below, the function has to be on.

2D Viewer

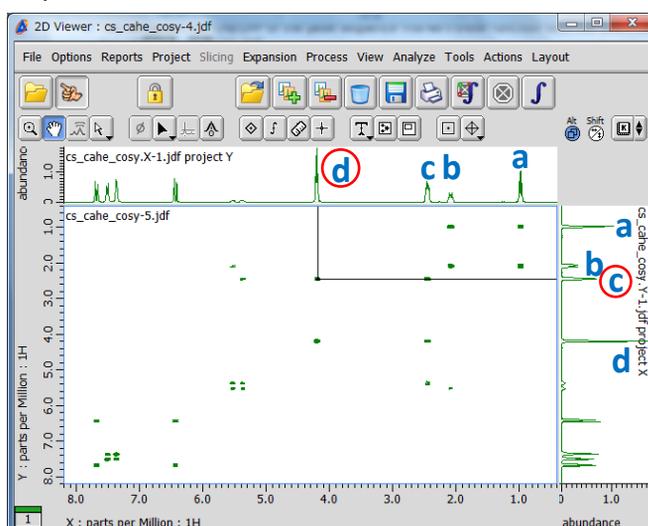



The peak sequence markers are shown as two perpendicular lines at **c** and **d** peaks

- 1 Push the **Peak** button  to activate the peak selection function.

Note that the cursor has changed into the **Peak symbol** .

- 2 Select a peak in the **2D Viewer** window with the cursor.

The peak sequence markers has been shown on the ^1H - ^1H COSY spectrum.

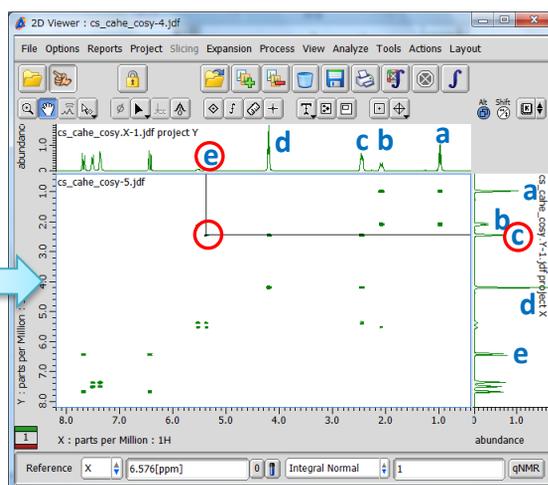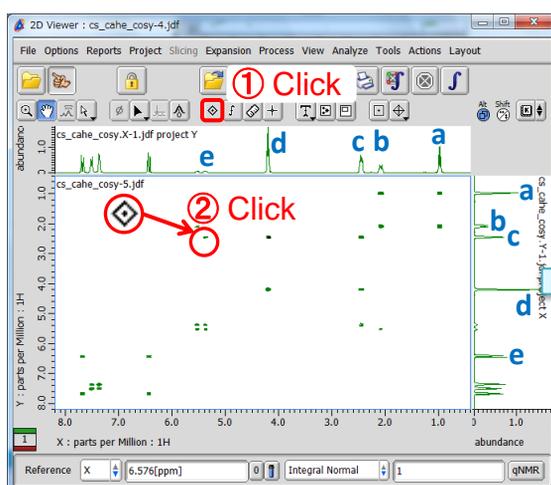


The peak sequence markers at **c** and **e** peaks

1/2

It is also possible to active/deactivate the peak sequence markers function as follows:

- ① Push and hold the right mouse button to display the context menu. Then select **Options — Peaks — Sequence Markers — Show Markers**.

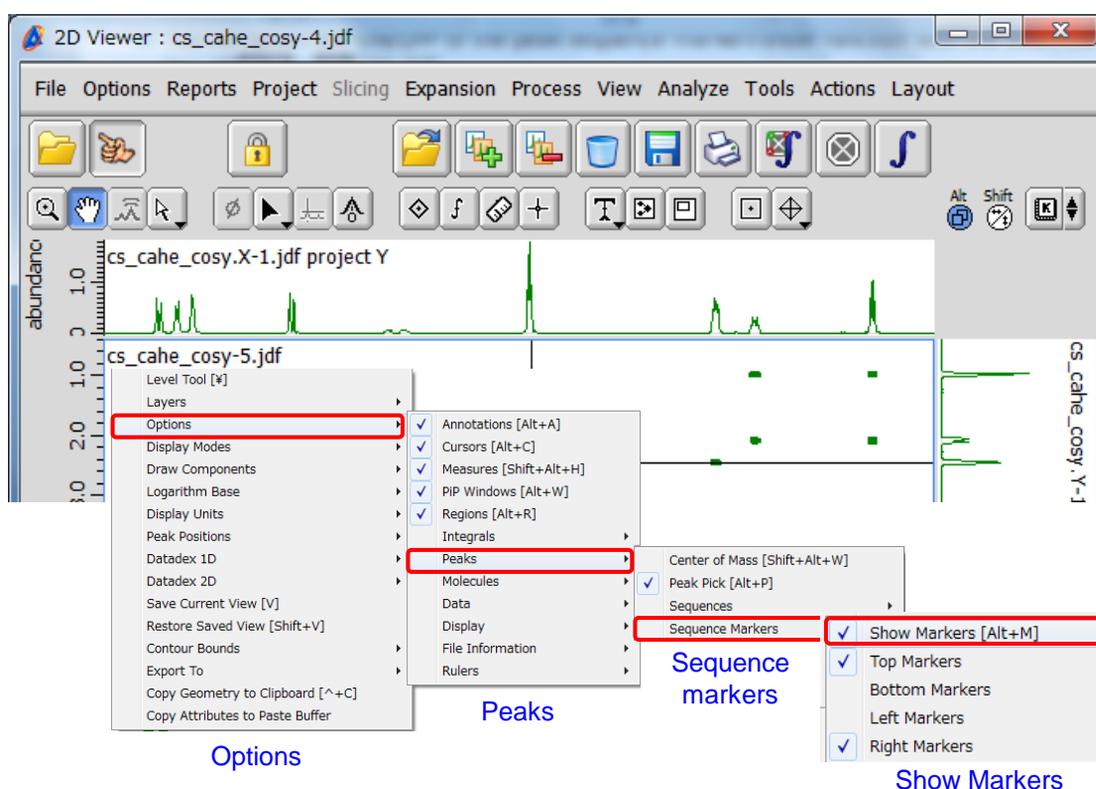

- ★ The keyboard shortcut (shortcut keys) is shown on the right.