

# Delta Tips

NMDT\_0052

## How to Delete Processed Data


NMR data processing software

Delta  
NMR Software  
v5.0



Delta software stores processed NMR data (i.e. Fourier transformed-data) in separate files apart from FID data. It is possible to quickly delete all files containing processed data while keeping the original FID data by using a function of the **File Browser** window.

- ★ Refer to User's Manual - Data Processing User's Manual for more detail on **File Browser** and file version numbers.

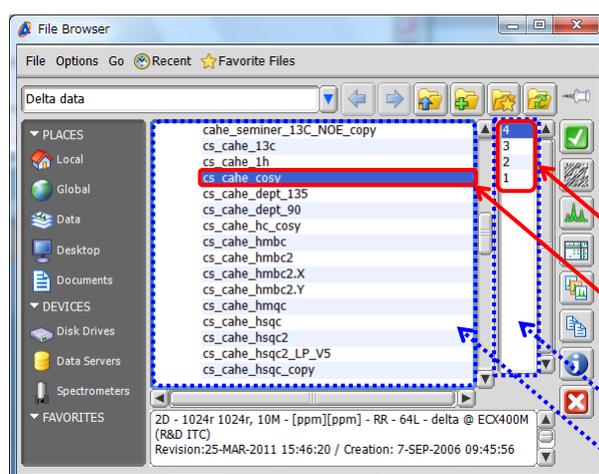

The **CS\_cahe\_COSY** data has been selected.

- ★ Delta typically stores FID data under version number 1 and FT-processed data under version numbers 2, 3, and so on. Each version number corresponds to one file.



The FID data (1) and the processed data (2-4) of the selected data file **cs\_cahe\_cosy**

Selected data file

Version Number Box

File List Box

Select a data file and go to **File – Purge Processed Versions** in the File Browser window. Note that all files with version numbers 2-4 have been deleted.

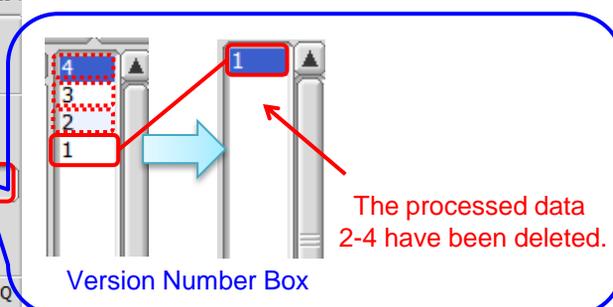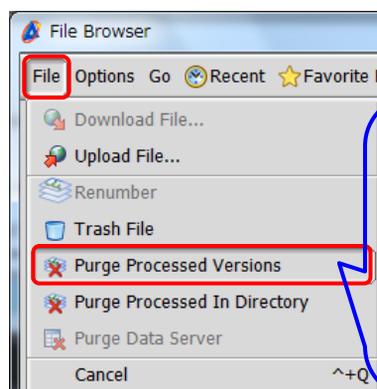

The processed data 2-4 have been deleted.

Version Number Box

- ★ The FID data (1) is still kept in the folder.
- ★ If you wish to delete one particular data file, select its version number in the version number box and select **File – Trash File** or push the **Delete** key.
- ★ If you wish to delete all processed data files in the folder, select **File – Purge Processed In Directory**.