

Delta Tips

NMDT_0054


Annotation Function

NMR data processing software

Delta
NMR Software
v5.0




Delta software has many functions and tools related to data processing and data presentation. In this issue, we will demonstrate the **Annotation** function.

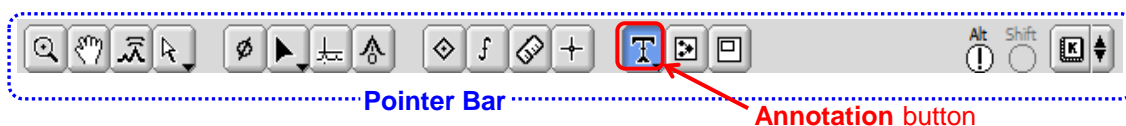




① Push the **Annotation** button  on the **Pointer Bar** in the Data Slate or Processor window.

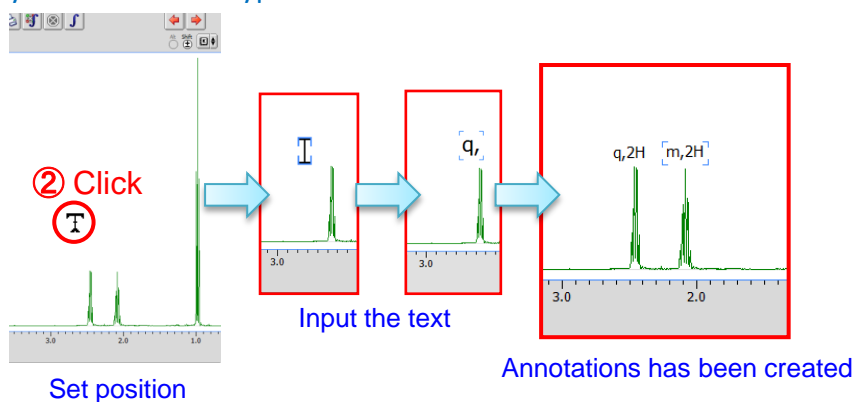

Note that the cursor has changed to the **Annotation** symbol .



② Select the place where you wish to create a text box with the cursor. An empty box has been displayed on the data. Type in the text of annotation.



★ You can display/hide annotations by pushing the  and  [Alt]+[A] keys.

It is also possible to create a **rectangle** or **arrow** on the data. Push and hold the **Annotation** button  to display the **Annotation** menu.

★ If you wish to create an arrow, select the **Arrow**  mode. Note that the cursor has changed into the **Arrow** symbol .

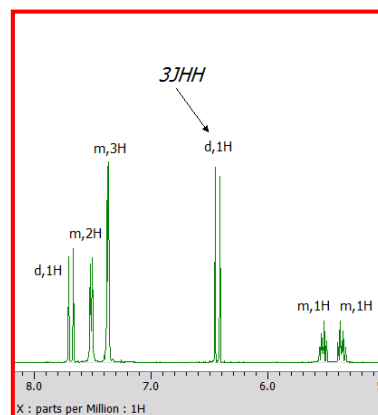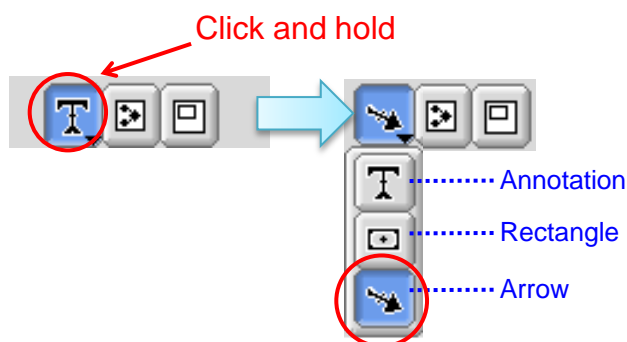

Annotations and arrow have been created