

Delta Tips

NMDT_0056

How to Customize Logo


NMR data processing software

Delta
NMR Software
v5.0



In Delta software, it is possible to replace the default JEOL logo by a logo (image) of your choice. In the example below, we will demonstrate how to change the logo by using the **Param Right** template and the **Page Layout Editor**.

★ For the details on **Page Layout Editor**, refer to NMDT_0055.

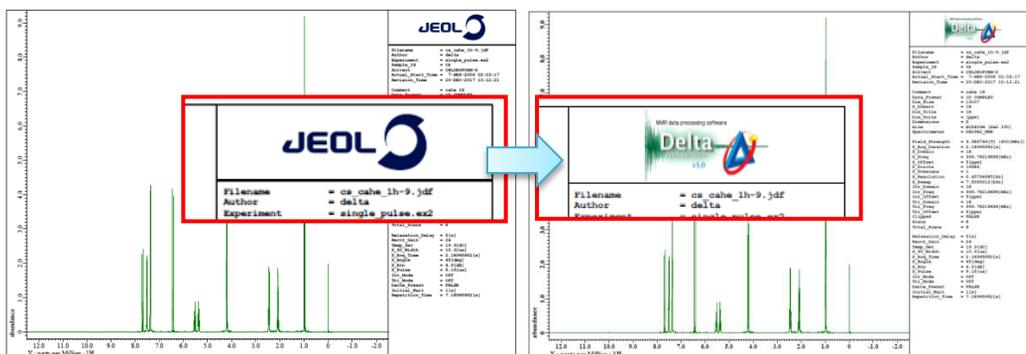

The default JEOL logo has been replaced by a Delta logo.

① Push the **Region** button  in the **Page Layout Editor** window.


Note that the cursor has changed into the **Region symbol** .

② Click the **Logo data area (Box 2)** with the cursor  to activate it.

③ Select **Import - EMF** in the **Page Layout Editor** window.

Select a logo file and push the **Open** button  in the **Specify emf file** window.

The selected logo has been loaded into **Box 2**.

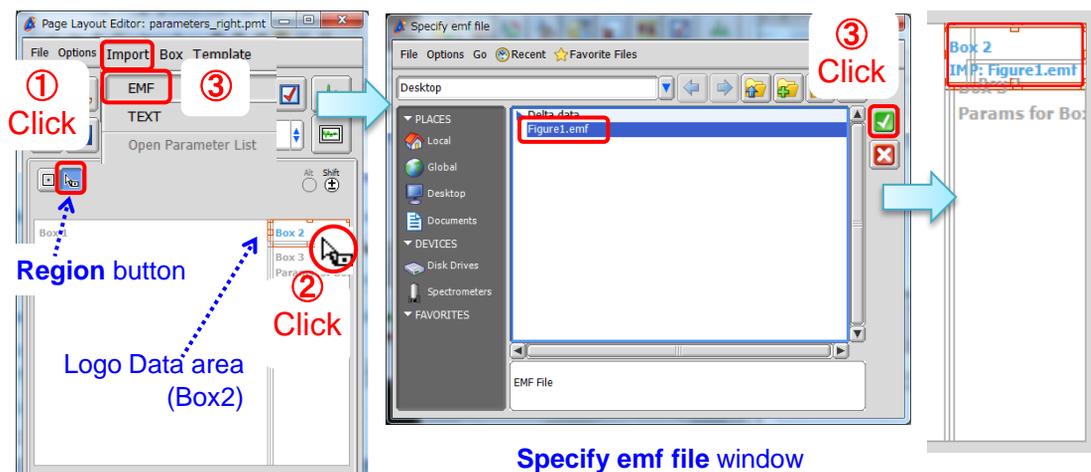

Specify emf file window


Figure1.emf file has been loaded into the Logo area.

1/2



④ Push the **Print Layout** button  to print the spectrum to a printer or PDF file.

Confirm that Figure1.emf is displayed in the logo area.

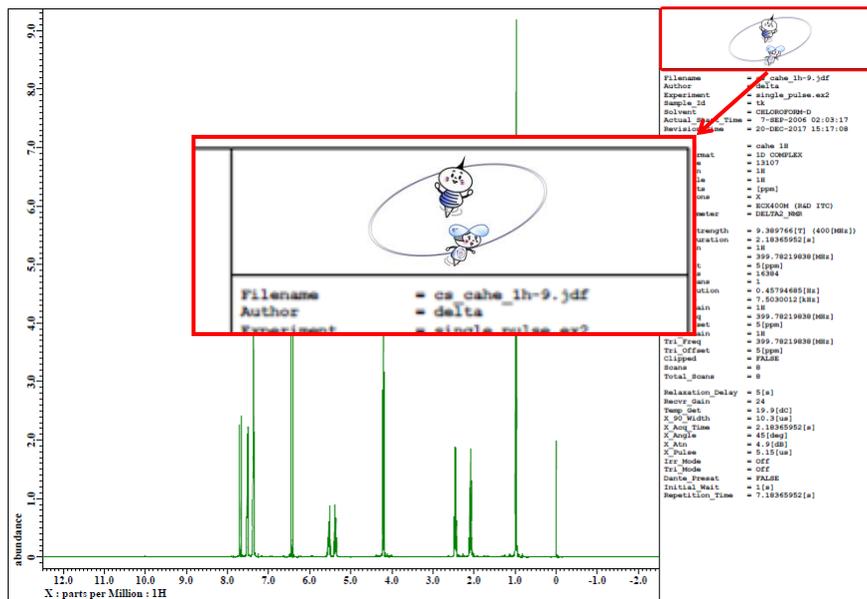

! Logo needs to be in the EMF (Enhanced MetaFile) format. If the logo of your choice is in another format, please convert it to EMF.

★ NMDT_0035 explains how to delete the logo.