

Delta Tips

NMDT_0069


How to Make a Print Layout (2)

NMR data processing software






< Customize lists of acquisition and processing parameters to print >

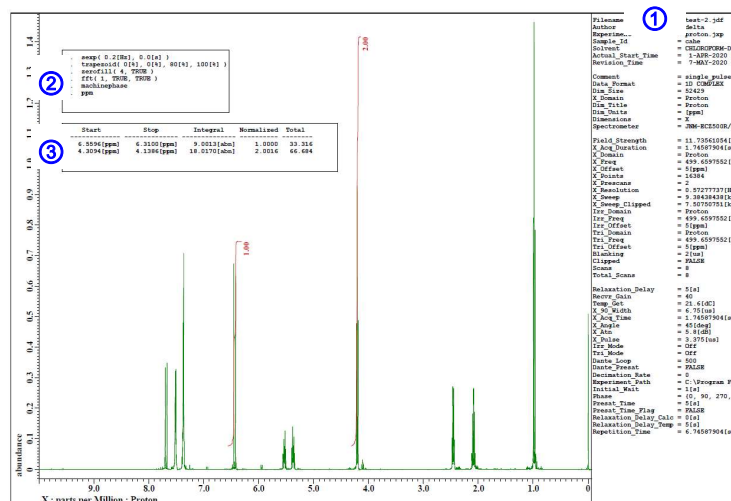In Delta software, it is possible to create sophisticated layouts for printing and presentation. This issue of Delta Tips explains additional functions of the **Page Layout Editor**. It is possible to customize various lists and text related to the sample and NMR data to print. In the examples below, we will customize lists of acquisition parameters ①, processing parameters ② and integration data ③.




★ For the details on creating various layouts, refer to the Delta Tips issue titled **How to Make a Print Layout (1)** (⇒NMDT_0068).


① Select a data. Select a data box and click the **Select Data** button  in the **Page Layout Editor** window. And then select the data with the cursor  to load this data to the data box. Note that the name of the selected data is displayed in the data box.

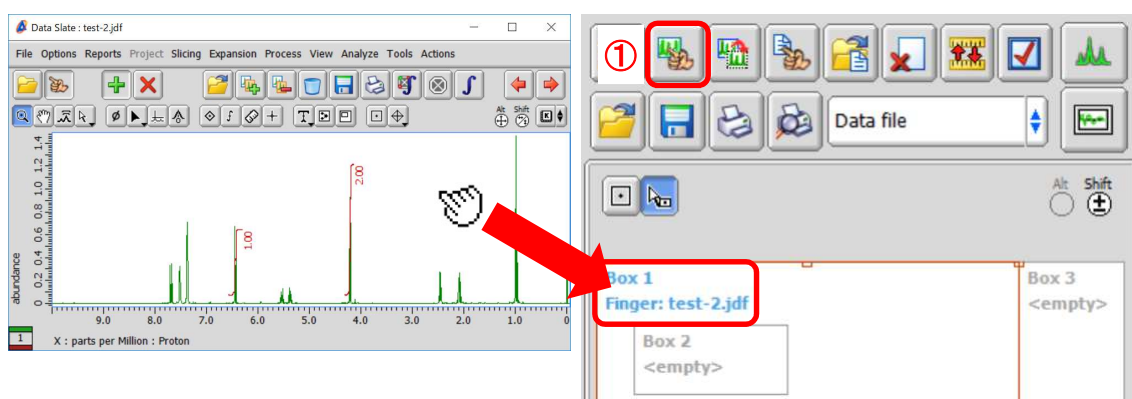


1/5

Delta Tips

- ★ If you wish to display lists of peaks and integrals, create the reports as follows:
Open a data and select **Reports—Peak Report** or **Integral Report** on the menu bar to create a peaks list or integrals list. Click the **File Information** button  in the **Delta Console** window to open the **File Information** window. Note that two reports have been shown in the **Reports** tab below. One of them reports peaks, the other one integrals.
- ★ For the details on creating and saving peak lists, refer to the Delta Tips issue titled **How to Create and Save Peak List** (⇒ NMDT_0014). For integral lists, refer to the Delta Tips issue titled **How to Create and Save Integral List** (⇒ NMDT_0015).

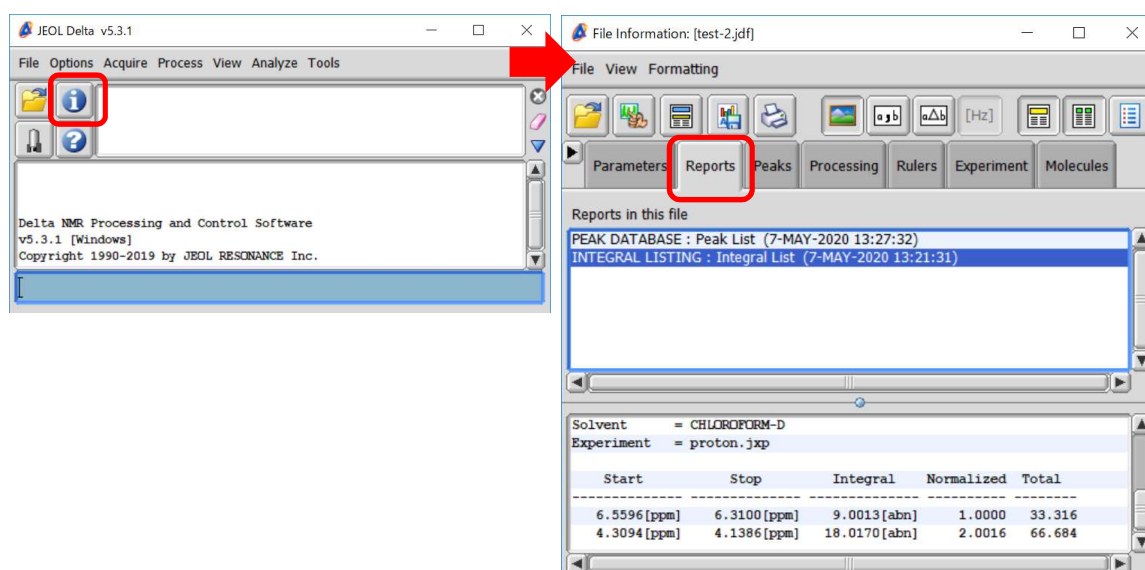

- ② Create a data box to load the report. Select the data box and select the source data in the list ② to link. In the example below, we have linked **Box 2** and **Box 3** to the source data **Box 1**.

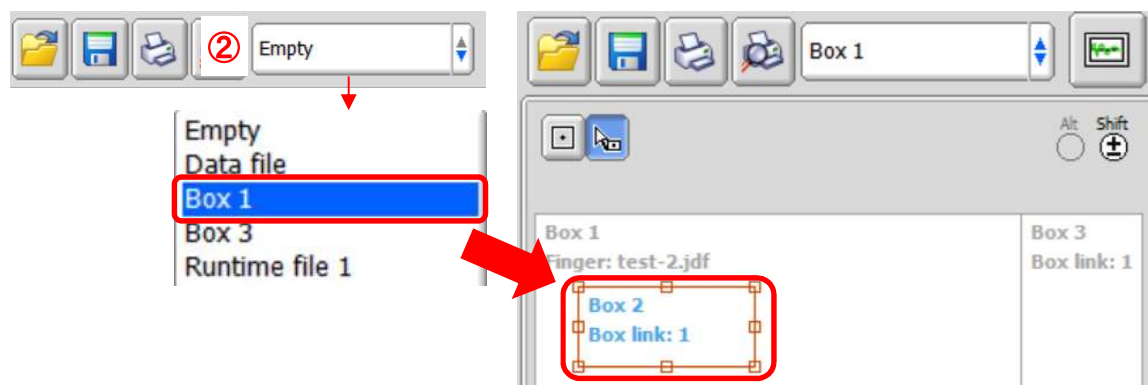


Delta Tips

NMR data processing software


Delta
NMR Software



- ③ Select the data box and click the button ③ to select **Data/ Parameters/ a Report/ a Processing List**. In the example below, we have set **Data** in **Box 1**, **Report** in **Box 2** and **Parameters** in **Box 3**.

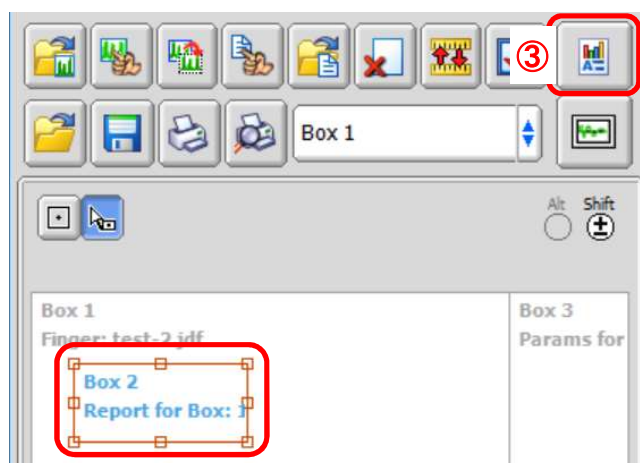

- ④ Select **Import—Select Report** to open the **Select Report** window. Select the report and click the **Select** button to load the report. In the example below, we have loaded an **Integral List** to **Box 2**.

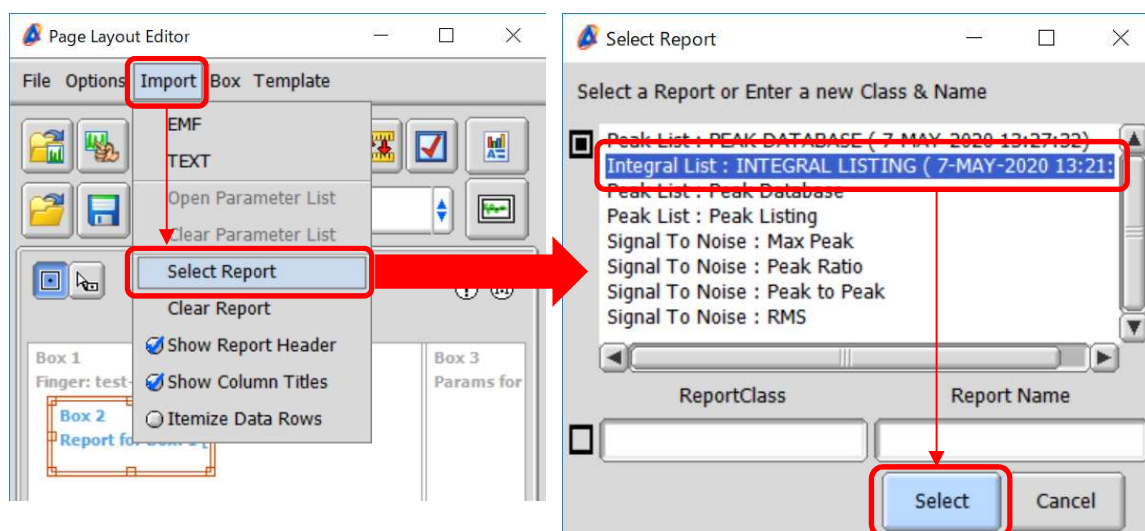




How to set font:


It is possible to set the size and style of letters in a data box as follows:
Select **Box — Attributes** to open the **Attributes** window.

- ★ For the details on the **Attributes** window, refer to the Delta Tips issue titled **Page Layout Editor (Part 1)** (⇒NMDT_0019).

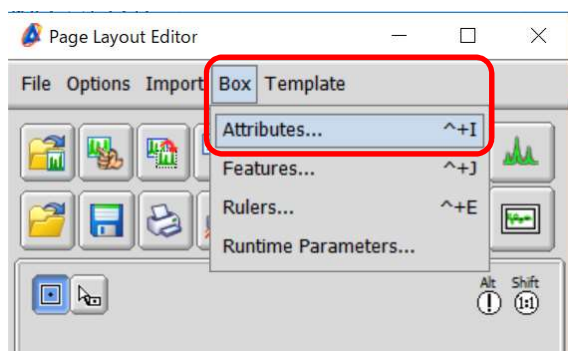

How to use templates:

It is possible to use preset templates as follows:

Select a template from the **Template** menu or click the **Open Template** button  to open the **Open File** window.

- ★ Refer to the Delta Tips issue titled **Page Layout Editor (Part 2)** (⇒NMDT_0055) for more details.

