

## **Cross Sections of Noodles Observed at Various Cooling Temperatures**

Instrument used: Scanning Electron Microscope (SEM)


SEM observation of a hydrated specimen requires the observation in a frozen state to preserve the morphology of the specimen. However, depending on the location of the SEM instrument, the use of liquid nitrogen may be difficult. In this article, cryo-techniques using various coolants were compared experimentally. The cross sections of noodles were observed using a low vacuum (LV) cooling holder under four different cooling conditions.

## **Comparison of cross sections of noodles at various temperatures**

Dried noodles were boiled under the same conditions. These noodles were then cooled at various temperatures, listed on the table below, before they were cut or fractured. The prepared specimens were observed by SEM in low vacuum (LV) mode.

Cooling temperature	(1) 4°C (in refrigerator)	(2) −20°C (in freezer)	(3) −79°C (with dry ice)	(4) –196°C (with liquid nitrogen)
Cross section preparation	Cross section fracture unit (CSC III)	Cross section fracture unit (CSC III)	Freeze fracture unit	Freeze fracture unit
Magnification x500 LV mode, uncoated LV cooling holder used				
Before observation	LV cooling holder was cooled in the refrigerator.	LV cooling holder was cooled in the freezer.	LV cooling holder was cooled with dry ice.	LV cooling holder was cooled with liquid nitrogen.
Status of cross section	Since the specimen was not frozen, a cut face was created by a razor blade. Due to the stickiness of the specimen, a clean cross section was not possible causing unclear boundaries between the starch and other parts.	Due to insufficient freezing of noodles, noodles were peeled away from the starch particles. The starch particle diameter was easy to measure.	Freeze fracturing was possible. Ice crystals around starch particles were confirmed before the increase of temperature.	Freeze fracturing was possible. When the freezing with liquid nitrogen was made, starch particles embedded in the ice crystals were confirmed.
Remarks	At 4°C, vaporization of water from the specimen occurred during SEM observation, allowing confirmation of the compositional distribution in the specimen. Detailed information on ice crystals was unable to be obtained.	Even at $-20^{\circ}$ C, vaporization of water from the specimen occurred during SEM observation. The cross section was non-uniform due to the various particle sizes of the starch.	At dry ice temperatures, the time to reach sublimation was very short. Observation in a frozen state was very difficult.	Stable observation was enabled with liquid nitrogen cooling.
Result	Impossible to fracture. The specimen was dehydrated during observation.	Impossible to freeze fracture. The specimen was dehydrated during observation.	Possible to freeze fracture. Observation in a frozen state was difficult. Observation in a freeze dried state was possible.	Possible to freeze-fracture. Observation in a frozen state was possible.

Conclusion: Liquid nitrogen is essential for SEM observation in a frozen state. Although the dry ice cooled specimen observation in a frozen state was difficult, necessary information can be obtained depending on some specimen type or observation purpose.

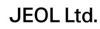


Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.

## Comparison of SEM images of noodles in frozen state and freeze dried state, using dry ice

In SEM observation using an LV cooling holder, the specimen temperature gradually increases and sublimation takes place, so the specimen state is transformed to a freeze dried state during the observation. In this state, specimen deformation can arise. In particular, at cooling with dry ice, the time to reach sublimation is very short and therefore, the specimen is likely to be observed in a freeze dried state.

In the frozen state, starch particles, embedded in ice crystals, are observed. In the freeze dried state, ice crystals around the starch particles are sublimated, revealing the emergence of voids.

In the freeze dried state, it is possible to confirm the size of starch particles, but difficult to view the status of the ice crystals.

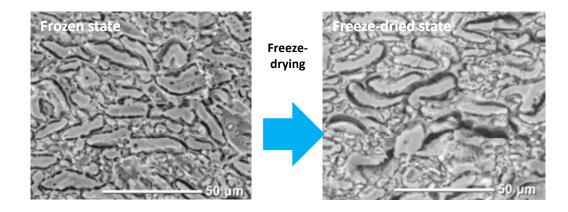

Copyright © 2017 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. LEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.



1-2 Musashino 3-chome Akishima Tokyo 196-8558 Japan Sales Division Tel. +81-3-6262-3560 Fax. +81-3-6262-3577 www.jeol.com ISO 9001 · ISO 14001 Certified

AUSTRALIA & NEW ZEALAND /JEOL(AUSTRALASIA) Pty.Ltd. Suite 1, L2 18 Aquatic Drive - Frenchs Forest NSW 2086 Australia • BELGIUM /JEOL (EUROPE) B.V. Planet II, Gebouw B Leuvensesteerweg 542, B-1990 Zaventem Belgium • BRAZIL /JEOL Brazil Instrumentos Cientificos Ltda. Av. Jabaquara, 2958 5<sup>s</sup> andar conjunto 52; 04046-500 Sao Paulo, SP Brazil • CANADA /JEOL CANADA, INC. 3275 1ere Rue, Local #8 St-Hubert, OC J3Y/8Y6, Canada • CHINA /JEOL(BELING) CO, LTD. Zhongkezyuan Building South Tower 27; Zhongguancun Nansanje Street No. 6, Haidian District, Beijing, PR.China • EGYPT /JEOL SERVICE BURRAU 3/dt Fl. Nile Center Bidg., Naval Street, Dokki, (Cairo), Egypt • FRANCE /JEOL (EUROPE) SAS Espace
Claude Monet, 1 Allee de Giverny 78290, Croissy-sur-Siene, France • GERMANY JGOL (GEHMANY) GMH Gui A Areger 30 8536 Freising, Germany • GRAET BITAIN & IRELAND /JEOL (UK.), LTD. JEOL House, Silver Court, Watchmead, Welwyn Garden City, Herts AL7 1LT, U.K. • ITALY /JEOL (ITALA) S.p.A. Palazzo Pacinotti - Milano 3 City, Va Ludovico II Moro, 6/A 20030 Basiglio(MI) Itay • KOREA /JEOL KOREA LTD. Dongwoo Bidg, 77; 1443, Prajae Daero, Gangdong-Gu, Seoul, 134-814, Korea • MALAVSIA) SIDU MALAVSIA) SIDU KALANSIA /JEOL (NALANSIA) SIDU (UK) A LUCIO: Caropration Road INS 77, Kelana Busina S 77, Kelana Java, 4730 Heating Java, Selangor, Malaysia • MEXICO /JEOL DE CV. Arkansas II Piso 2 Colonia Napoles Delegacion Benito Juarez, C.P. 03810 Mexico J FL, LTD. 2 Coropration Road IIO-12 Corporation Place SingApper 6 18494 • TAIWAN /JEOL (CHONC P) LEOL (Scandinaviska)AB Hammarbacken 6A, Box 716, Bid Z. 127473, Moscow, Russian Federation • SCANDINAVIA /SWEDEN JEOL (Skandinaviska)AB Hammarbacken 6A, Box 716, Bid Z. J127473, Moscow, Russian Federation • SCANDINAVIA /SWEDEN JEOL (Skandinaviska)AB Hammarbacken 6A, Box 716, Bid Z. J127473, Moscow, Russian Federation • SCANDINAVIA /SWEDEN JEOL (Skandinaviska)AB Hammarbacken 6A, Box 716, Bid Z. J127473, Moscow, Russian Federation • SCANDINAVIA /SWEDEN JEOL (Skandinaviska)AB Hammarbacken