

Application of ROYALPROBE™ P+ to catalyst

Product used : Nuclear Magnetic Resonance (NMR)

Feature of ROYALPROBE™ P+

The ROYALPROBE™ P+ is a newly developed solution NMR triple-resonance probe that is capable of ^1H , X, and ^{31}P triple-resonance measurements. It can be used as a triple-resonance probe even on a standard 2-channel ECZ Luminous™ console due to the MFDS function¹⁾. Fig. 1 is a wiring diagram of the ROYALPROBE™ P+ connected to a 2-channel ECZ Luminous™ console. In the case of conventional triple-resonance probe, it is necessary to manually exchange the corresponding frequency filter every time the X nucleus is changed, therefore it is not possible to change the X nucleus by AUTO TUNING UNIT only. In contrast, the ROYALPROBE™ P+ uses a newly developed PX diplexer that can serve as a filter for various X nuclei. This allows us to tune the probe on all three channels automatically using the AUTO TUNING UNIT. Furthermore, compared to a conventional triple-resonance probe, the sensitivity of ^{13}C , ^{31}P and other nuclei has been greatly improved, making it much more versatile. Here we demonstrate an application of this new probe to a sample of catalyst containing ^{31}P .

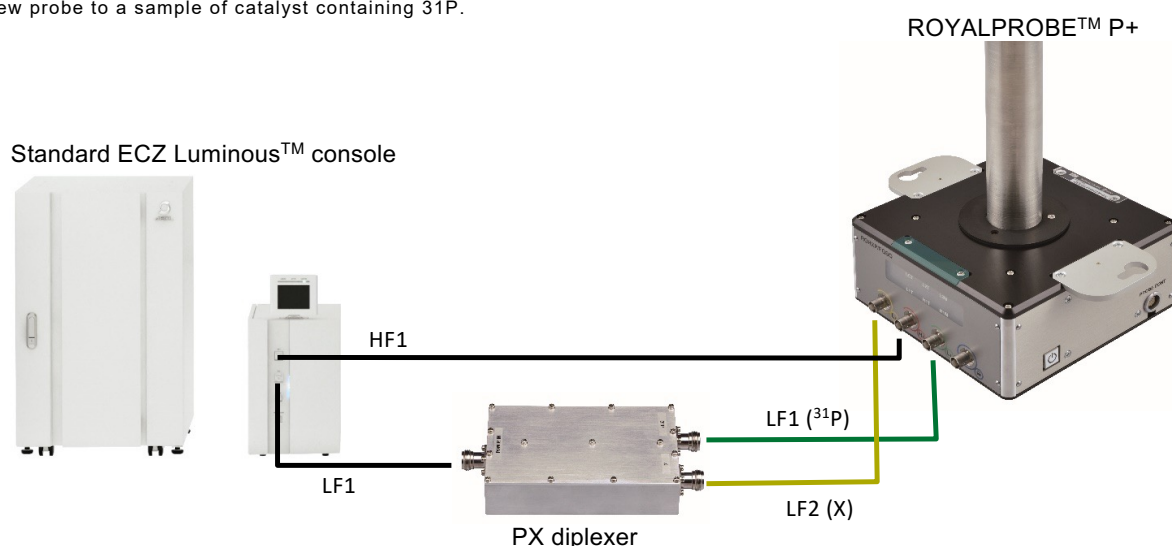


Fig. 1: Wiring of ROYALPROBE™ P+ and PX diplexer when using ECZ Luminous™ in standard configuration

Triple-resonance experiment of ^1H , ^{11}B and ^{31}P

For many organic compounds, the ^1H chemical shift range is about 10 ppm, but it is known that the ^1H chemical shift may be significantly shifted by metals. Fig. 2 shows a ^1H NMR spectrum of Ru-MACHO®-BH* which is known as an ester hydrogenation catalyst. A very characteristic ^1H signal can be observed in high field above -10 ppm which is a very unusual value. In addition, a very broad and split signal at -2.3 ppm is observed. Since this sample contains heteroatoms such as ^{31}P and ^{11}B that affect the shape of ^1H signal, decoupling of these nuclides can help us identify the ^1H nuclei close to the phosphorus and boron atoms.

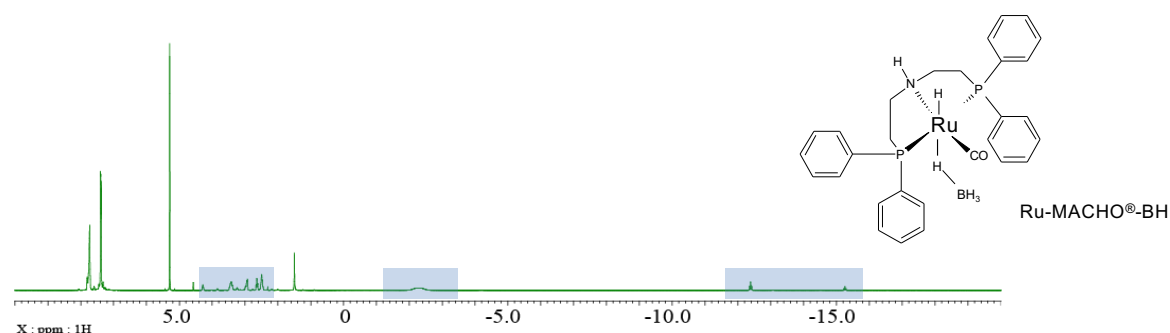


Fig. 2: ^1H NMR spectrum of 3mg Ru-MACHO®-BH in CD_2Cl_2

*Ru-MACHO® is a registered trademark of TAKASAGO INTERNATIONAL CORPORATION.

The spectrum shown in Fig. 2 is further expanded in Fig. 3-5. The CH₂ range is shown in Fig. 3. The signal of BH₄ group is shown in Fig. 4. Finally, the signals of Ru-H protons are shown in Fig. 5. a) refers to standard ¹H spectra, b) refers to ¹H{¹¹B} spectra, c) refers to ¹H{³¹P} spectra, and d) refers to ¹H{¹¹B}{³¹P} spectra. By careful comparison of the coupled and uncoupled ¹H spectra, it is possible to identify the signals of the CH₂ and Ru-H groups and the signal of the BH₄ group. In catalysts containing transition metals, such as this sample, it is common for various stereoisomers to exist and be present. For this reason, more signals are often observed than one would expect from the structural formula. This phenomenon can also be confirmed in Fig. 3-5 where two sets of signals are clearly observed. Therefore, narrowing and simplifying signals by heteronuclear decoupling may help reduce overlaps and identify minor components and/or impurities.

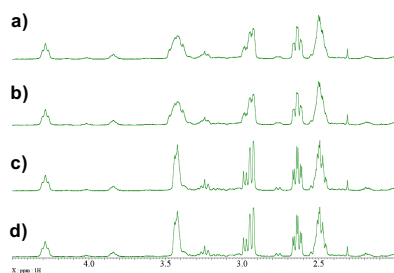


Fig. 3: Expanded region of CH₂ groups

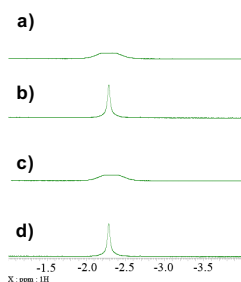


Fig. 4: Expanded region of BH₄ group

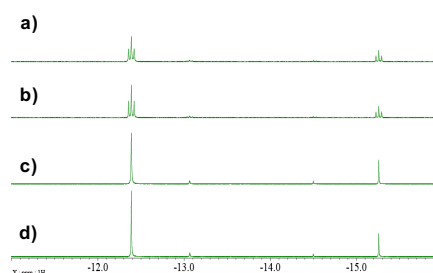


Fig. 5: Expanded region of Ru-H group

Triple-resonance experiment of ¹³C, ¹H and ³¹P

Fig. 6 shows a standard ¹³C{¹H} spectrum of the catalyst, whilst Fig. 7 shows the expansion indicated by the blue rectangle in Fig. 6. As you can see in the ¹³C{¹H}{³¹P} spectrum (Fig 7b), the signals of CH groups and quaternary carbons in the vicinity of ³¹P are very easy to distinguish by comparing the double-resonance and triple-resonance spectra. As ¹³C is a low sensitivity nuclide due to its low natural abundance (1.1%), the signals of carbon atoms coupled with ³¹P may be difficult to detect in the case of diluted samples like this one. For this reason, ³¹P-decoupling simplifies the spectrum and allows us to observe signals which could be probably missed.

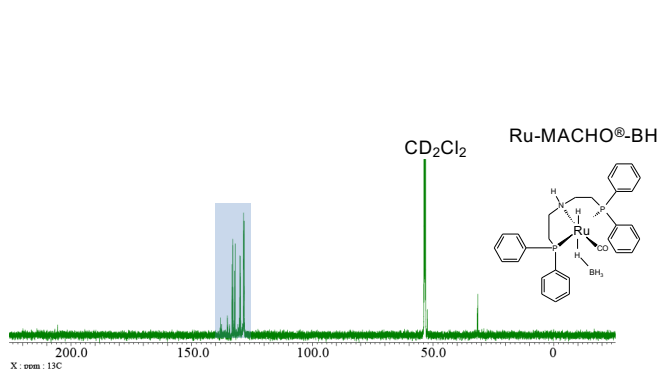


Fig. 6: ¹³C{¹H} spectrum

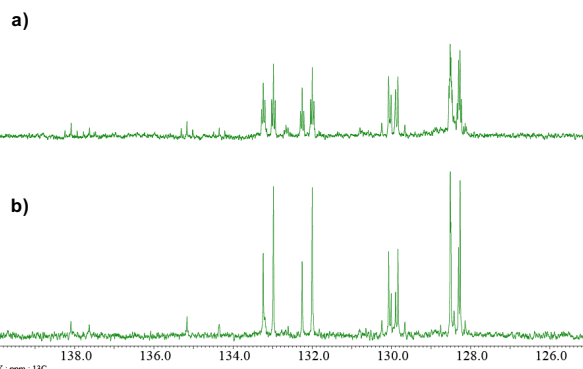


Fig. 7: Expanded aromatic region of a) ¹³C{¹H}, b) ¹³C{¹H}{³¹P} spectra

In the case of samples containing stereoisomers or multiple components, complex spectral patterns are often observed. Therefore, the capability to decouple ³¹P and other heteronuclei and to employ advanced correlation experiments is very important in the analysis of complex samples. Therefore, the ROYALPROBE™ P+ and other triple-resonance probes utilizing the MFDS functionality of the ECZ Luminous™ are of great help to those who synthesize and analyze such samples.

Reference

1) Application Note NM220004E

Sample courtesy of TAKASAGO INTERNATIONAL CORPORATION



3-1-2 Musashino Akishima Tokyo 196-8558 Japan Sales Division Tel. +81-3-6262-3560 Fax. +81-3-6262-3577
www.jeol.com ISO 9001 • ISO 14001 Certified

• AUSTRALIA & NEW ZEALAND • BELGIUM • BRAZIL • CANADA • CHINA • EGYPT • FRANCE • GERMANY • GREAT BRITAIN & IRELAND • ITALY • KOREA • MALAYSIA • MEXICO • RUSSIA • SCANDINAVIA
• SINGAPORE • TAIWAN • THE NETHERLANDS • USA



Copyright © 2022 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.